



GDM-49-SCI-HC~TSL01



T-49-SCI-HC

⚠ WARNING!



Be sure to read and fully understand this document before installing, operating, maintaining or servicing this appliance. Failure to do so can result in appliance failure, property damage, serious injury or death. Appliance failure, injury or property damage due to improper installation is not covered by warranty.

📌 USER ACTION!

TRUE tracks the history of your appliance by its serial number. For easy reference, record your appliances full model name and serial number below. This information is on the serial label. Serial label location varies by appliance.

Model Name:

Serial Number:

True Manufacturing Co., Inc.

2001 East Terra Lane • O'Fallon, Missouri 63366-4434
(636) 240-2400 • FAX: (636)-272-2408

International FAX: (636)-272-7546 • (800)-325-6152

Parts Department: (800)-424-TRUE (424-8783)
Parts Department FAX: (636)-272-9471



INSTALLATION MANUAL

GDM/T-Series Scientific

Original Instructions

North America – Canada and Caribbean

Warranty Phone +1 855-878-9277
Warranty Fax +1 636-980-8510
Warranty Email warrantyinquiries@truemfg.com
Technical Phone +1 855-372-1368
Technical Email service@truemfg.com
7:00 am–6:00 pm CST Monday–Friday,
8:00 am–12:00 pm Saturday

Mexico

Phone +52 555-804-6343/44
service-mexicocity@truemfg.com
9:00 am–5:30 pm M–F

Latin America

Phone: +52 555-804-6343/44
servicelatam@truemfg.com
9:00 am–5:30 pm M–F

UK, Ireland, Middle East, Africa & India

Phone: +44 (0) 800-783-2049
Service-emea@truemfg.com
8:30 am–5:00 pm M–F

Australia

Phone: +61 2-9618-9999
service-aus@truemfg.com
8:30 am–5:00 pm M–F

European Union & Commonwealth of Independent States

Phone: + 41 61 563 0705
service-emea@truemfg.com
8:00 am–5:00 pm M–F



THANK YOU

FOR YOUR PURCHASE

Congratulations!

The primary purpose of this document is to assist the installation, maintenance, and servicing of your TRUE appliance. This document contains information important to safety, operation, maintenance, and servicing. **DO NOT** discard this document. TRUE is solely the appliance manufacturer. For assistance locating a refrigeration service technician in your area for installation, servicing or maintenance, please visit our Service Company Locator at truemfg.com/support/service-locator.

NOTICE!

Your appliance may not exactly match the figures shown in this manual.

Contents






Preface	
Signal & Symbol Definitions	3
Important Safety Information	
Basic Safety & Operation Warnings	4
Personal Injury Warnings	5
Appliance Disposal Warnings	6
Hydrocarbon Refrigerant Warnings	6
Electrical Safety Warnings	7
About Your Appliance & Installation Requirements	
Appliance Specifications	8
Clearances	9
Electrical Requirements	10
Electrical Installation & Safety	10
Grounded Electrical Duplex with Cover	11
Installation & Setup	
Uncrating	12
Appliance Location	15
Leveling Screw, Leg or Castor Installation	15
Level the Appliance	17
Seal the Appliance to the Floor	17
Shelf Installation	18
Chromatography Accessories	19
Appliance Operation	
Startup	20
Temperature Control & Light Switch Location	21
General Sequence of Operation	23
Temperature Control	25
Electronic Temperature Control	
Lock / Unlock the Control	27
Turn Off / On the Control	27
Change the Set Point	28
Initiate Manual Defrost	28
Change System of Measure	29
Display Probe Temperature	30
Adjust Temperature Alarm Ranges	31
Unlock the Control	33
Turn Off the Control	34
Turn On the Control	34
Change the Set Point	35
Initiate Manual Defrost	35
Change System of Measure	36
Display Probe Temperatures	36
Temperature Alarms	37
Door Open Alarm	38
Adjust Temperature Alarm Ranges	38
High (thi) & Low (tLo) Temperature Log	41
Maintenance & Servicing	
Component Replacement	43
Recommended Maintenance	44
Condenser Coil Cleaning	46
General Surface Care & Cleaning	48
Stainless Steel Care & Cleaning	49
8 Tips to Help Prevent Rust on Stainless Steel	50
Slide Door Removal	51
Slide Door Adjustment	53
Slide Door Hold-Open Feature	53
Warranty	
Warranty Information	54











Preface

The warning, guidelines, and recommendations within this document are meant to prevent appliance damage, property damage, injury, or death. Please carefully read all warnings, guidelines, and recommendations before proceeding to ensure the continued safe use and maintenance of your TRUE appliance.

Signal & Symbol Definitions

Below are symbols you may see in this document. Some symbols may not appear.

Signal Word Definitions	
 DANGER!	An imminently hazardous situation which, if not avoided, will result in serious injury or death.
 WARNING!	A potentially hazardous situation which, if not avoided, can result in serious injury or death.
 CAUTION!	A potentially hazardous situation which, if not avoided, may result in minor or moderate injury; an unsafe practice.
 USER ACTION!	User action alert, follow all recommendations to avoid appliance or product damage.
 NOTICE!	Important information not related to hazards or risk of personal injury.

Safety Symbols	
	Safety alert; alerts reader to potential physical injury hazards. Obey all safety messages following this symbol to avoid possible injury or death.
 	Flammable material; fire or explosion hazard.
	Electrical shock hazard.
	Tipping hazard; tip-over hazard.
	Sharp element; cut or sever hazard.
	Eye hazard; risk of eye injury.
	Exploding pressurized cylinder hazard.
	Slippery surface hazard.
	Crush or cut hazard.

Additional Symbols	
	Mandatory action alert symbol; alerts reader to required or recommended actions. Obey all messages and recommendations following this symbol to avoid appliance or product damage.
	Important information not related to hazards or risk of personal injury.
	Review and understand the installation manual before installing, operating, or servicing.
	Wear eye protection.
	Wear protective gloves.
	Secure gas cylinders to prevent falling cylinders.
	DO NOT use extension cord.
	DO NOT use adaptor plugs.
	DO NOT dispose of with other household waste.



Important Safety Information

Important Safety Information

Basic Safety & Operation Warnings

Follow basic safety precautions, including the following, to reduce risk of personal injury, electric shock, fire, or death.


⚠ WARNING!

 	<p>Be sure to read and fully understand this document before installing, operating, maintaining, or servicing this appliance. Failure to do so can result in appliance failure, property damage, serious injury, or death. Appliance failure, personal injury, or property damage due to improper installation is not covered by warranty.</p> <ul style="list-style-type: none"> Failure to install, operate, and maintain the appliance as detailed in this document will negatively affect safety, appliance performance, component life, and warranty coverage. Only qualified technicians should install and service the appliance. For assistance locating a refrigeration service technician in your area for installation, servicing or maintenance, please visit our Service Company Locator at truemfg.com/support/service-locator. TRUE is solely the appliance manufacturer and is not responsible for installation. This appliance is not to be used, cleaned, or maintained by persons (including children) with reduced physical, sensory, or mental capabilities or lack of experience and knowledge, without proper supervision or instruction concerning the appliance by a person responsible for their safety. DO NOT install or operate equipment that has been misused, abused, neglected, damaged, or altered/modified from original manufactured specifications. DO NOT modify or alter the appliance. DO NOT use electrical appliances inside the food storage compartments of the appliance unless the appliances are approved by the manufacturer. The appliance owner is responsible for performing a Personal Protective Equipment (PPE) Hazard Assessment and to ensure adequate protection during maintenance and cleaning procedures. Use appropriate tools, safety equipment, and PPE during installation and servicing. Only use the appliance for its intended purpose as described in this document. All shelving must be properly installed and adhere to load limits. Incorrect installation and improperly loaded or overloaded shelves can result in appliance damage, product damage, or personal injury. Keep the area surrounding the appliance clean and dry to avoid personal injury or appliance damage from debris or pests.
--	---

ⓘ USER ACTION!

	<ul style="list-style-type: none"> The appliance must be installed in accordance with all applicable laws, codes, and regulations. This appliance is to be installed in accordance with the Safety Standard for Refrigeration Systems, ANSI/ASHRAE 15.
--	--

ⓘ NOTICE!

	<p>The manufacturer is not responsible for injury or damage resulting from improper, incorrect, and unreasonable use.</p>
---	---

Important Safety Information (cont.)

Personal Injury Warnings

⚠ DANGER!





	DO NOT allow children to play with or in the appliance. Child entrapment or personal injury can occur.
	<p>DO NOT store or use the following in the vicinity of this or any other appliance:</p> <ul style="list-style-type: none"> Gasoline or other flammable vapors and liquids Combustible or explosive substances, such as aerosol cans with a flammable propellant Other volatile or flammable substances
 	<p>Contact TRUE Manufacturing to locate refrigerant lines and electrical wiring before drilling, cutting or puncturing interior or exterior walls. Failure to do so could result in damage, personal injury, or death.</p>

⚠ WARNING!

	<p>Only qualified technicians should install and service the appliance. For assistance locating a refrigeration service technician in your area for installation, servicing or maintenance, please visit our Service Company Locator at truemfg.com/support/service-locator. TRUE is solely the appliance manufacturer and is not responsible for installation.</p> <ul style="list-style-type: none"> Use appropriate tools, safety equipment, and personal protective equipment (PPE) during installation and servicing. DO NOT touch the cold surfaces in the freezer compartment when hands are damp or wet. Skin may stick to extremely cold surfaces.
	<p>This product can expose you to chemicals including Di-(2-Ethyl hexyl) Phthalate (DEHP), which is known to the State of California to cause cancer and birth defects or other reproductive harm. For more information go to P65warnings.ca.gov.</p>
	<p>Slippery Surfaces! Moisture from improper drainage can create slippery surfaces near the appliance. It is your duty to immediately warn your customers of, and dry, the slippery surface. All wet floor areas must be marked with a wet floor sign.</p>
	<p>Sharp edges! Take care when moving, installing, cleaning, servicing, and maintaining the appliance to avoid cuts. Be sure to take care when reaching under the appliance or handling metal components.</p> <ul style="list-style-type: none"> Keep fingers out of pinch point areas, such as the space between appliance doors and surrounding cabinetry. Take care closing doors with children nearby.
	<p>Tip over hazard! Appliance may pose a tipping hazard when uncrating, installing, or moving the appliance. Take appropriate safety precautions. Use of tip over restraints may only reduce (not eliminate) the tipping hazard. Never allow children to climb or hang on drawers, doors, or shelves.</p>
	<p>Crush or cut hazard! Keep clear when uncrating, installing, moving, or servicing the appliance.</p>
	<p>Risk of electric shock or burn! See "Electrical Safety Warnings" for more information.</p>


Important Safety Information (cont.)

Appliance Disposal Warnings

⚠ DANGER!	
	<p>Risk of child entrapment!</p> <p>Children can get trapped inside discarded appliances and suffocate. Never dispose of your appliance without taking precautions to prevent child entrapment, even if the appliance only sits unattended for a short period of time.</p> <p>Child entrapment precautions include the following:</p> <ul style="list-style-type: none"> • Remove all doors (or drawers for drawer appliances). • Leave all interior drawers and shelving in place to make climbing (and fitting) inside the appliance more difficult.
	
	<p>Risk of fire or explosion! Flammable refrigerant and insulation used. Dispose of the appliance in accordance with all applicable laws, codes, and regulations. Follow all safety precautions associated with handling flammable refrigerant and insulation.</p>
	<p>DO NOT dispose of your appliance with household waste.</p>

Hydrocarbon Refrigerant Warnings







TRUE appliances use hydrocarbon refrigerant (R-290/513A/600a). Check the serial label to identify the appliance's refrigerant. Serial label location varies by model.

⚠ DANGER!	
	<p>Risk of fire or explosion! Flammable refrigerant used.</p> <ul style="list-style-type: none"> • All servicing and maintenance must be performed by qualified technicians to minimize the risk of fire or personal injury due to incorrect parts or improper service. • Check the serial label to identify the appliance's refrigerant. Serial label location varies by model. • DO NOT damage the refrigeration system during transportation and installation. • If the appliance is damaged, check the refrigeration system for damage before installing or putting into service. • Never use sharp objects or tools to remove ice or frost. DO NOT use mechanical devices to accelerate defrost. • Dispose of the appliance in accordance with all applicable laws, codes, and regulations. Follow all safety precautions associated with handling flammable refrigerant. • DO NOT use electrical appliances inside the food storage compartments of the appliances unless the appliances are of the type recommended by the manufacturer.


Important Safety Information (cont.)

Electrical Safety Warnings

⚠ DANGER!

     	<p>Risk of electric shock, burn, or fire! Failure to comply with these electrical warnings may result in appliance damage, property damage, electrical shock, burns, or fire leading to serious personal injury or death.</p> <ul style="list-style-type: none"> It is the appliance owner's responsibility to ensure the electrical connection meets all applicable building codes. Before connecting the appliance to the power supply, verify the supply voltage and circuit rating match the appliance. Correct improper supply voltage or circuit size immediately. Before connecting the appliance to the power supply, verify the power supply is correctly grounded. If the power supply is not grounded, correct immediately. TRUE recommends hiring a qualified electrician to inspect the wall outlet and circuit to ensure they are properly grounded. DO NOT clean appliance with a pressure washer or hose. DO NOT immerse power cord in water. Always unplug the appliance or disconnect power before installation or servicing. Powering off an electronic control or setting temperature controls to 0 (off position) does not remove power from all components. The appliance should receive power from its own individual dedicated electrical circuit to avoid overloading the power supply. The original equipment manufacturer (OEM) power cord has a grounding plug to minimize the possibility of electrical shock. <ul style="list-style-type: none"> Never remove the ground prong from the power cord! For personal safety, this appliance must be properly grounded. Never use an extension cord! An extension cord is any component that adds length to the OEM power cord when connecting the cord to a power source. Never use an adapter plug! An adapter plug alters the OEM plug configuration when connecting the plug to a power source. Never use a power cord that shows cracks or abrasion damage along its length or at either end. <ul style="list-style-type: none"> Immediately have a qualified technician replace damaged OEM power cords with OEM components. Never unplug the appliance by pulling on the power cord! Always grip the plug and pull the plug straight out of the outlet. Take care to prevent rolling over or damaging the power cord when moving the appliance. Never use the power cord to prevent appliance movement! Always use adequate means to keep the appliance in place without transmitting stress to the power cord. Keep power cord away from heated surfaces. DO NOT let power cord hang over edge of table or counter. DO NOT excessively bend or place heavy articles on the power cord.
--	---

ⓘ NOTICE!

	<p>TRUE will not warranty the following:</p> <ul style="list-style-type: none"> Compressor failures due to improper incoming voltage. Appliance with tampered OEM power cords. Appliance connected to adapters or extension cords. <p>For more details, see TRUE's full warranty statement.</p>
---	---

About Your Appliance & Installation Requirements

About Your Appliance & Installation Requirements

! USER ACTION!



TRUE is not responsible for damage incurred during shipment. Always carefully inspect for freight damage before receiving and installing the appliance. If there is damage, note all damage on the delivery receipt, immediately file a claim with the delivery freight carrier, and contact TRUE. **Do not install the appliance or put it in service.**

Thank you for choosing TRUE Manufacturing to meet your refrigeration needs. TRUE highly recommends a qualified technician and electrician install your appliance to ensure correct installation. The cost of professional installation is money well spent. Only qualified technicians should install and service the appliance.

For assistance locating a refrigeration service technician in your area for installation, servicing or maintenance, please visit our Service Company Locator at truemfg.com/support/service-locator. TRUE is solely the appliance manufacturer and is not responsible for installation.

The appliance owner is responsible for proper installation and maintaining the appliance as described in this document. Routine care and maintenance procedures are not covered by TRUE's warranty.

Appliance Specifications

Some things to know about your appliance are as follows:

- Appliance is not for the storage and/or display of potentially hazardous foods when the temperature control is set above 41°F (5°C).
- Appliance is not suitable for an area where a pressure washer or hose may be used.
- Always plug the appliance into its own individual dedicated electrical circuit!
- **DO NOT** use extension cords or adapter plugs.
- Before connecting your appliance to the power supply, verify the incoming voltage ($\pm 5\%$) and the amps match the operation ratings on the appliance's serial label. Correct improper incoming voltage or amps immediately. Serial label location varies by model.
- Before connecting your appliance to the power supply, verify the power supply is correctly grounded. If the power supply is not grounded, correct immediately.
- Ensure the installation location will provide adequate clearances and sufficient airflow for the appliance. See "Clearances" (pg. 9).
- Read and follow all warnings and maintenance instructions. Failure to do so may result in damage and void the warranty on your appliance.

About Your Appliance & Installation Requirements (cont.)

Clearances

! NOTICE!	
!	<p>Warranty is void if ventilation is insufficient.</p> <p>Be sure your appliance has the required surrounding clearances for ventilation purposes. Keep all ventilation openings in the appliance enclosure or structure housing the appliance clear of obstruction.</p>

MODEL	TOP	SIDES	BACK
GDM Freezer	3" (76.2 mm)	0" (0 mm)	3" (76.2 mm)
GDM Refrigerator	0" (0 mm)	0" (0 mm)	1" (25.4 mm)
T-Series Freezer	3" (76.2 mm)	0" (0 mm)	3" (76.2 mm)
T-Series Refrigerator	0" (0 mm)	0" (0 mm)	1" (25.4 mm)

About Your Appliance & Installation Requirements (cont.)

Electrical Requirements

ⓘ USER ACTION!




 Find a copy of the wiring diagram with our serial number lookup at truemfg.com/support/serial-number-lookup.

Wire gauge chart (115V)												
115 Volts	Distance in Feet to Center of Load											
AMPS	20	30	40	50	60	70	80	90	100	120	140	160
2	14	14	14	14	14	14	14	14	14	14	14	14
3	14	14	14	14	14	14	14	14	14	14	14	12
4	14	14	14	14	14	14	14	14	14	12	12	12
5	14	14	14	14	14	14	14	12	12	12	10	10
6	14	14	14	14	14	14	12	12	12	10	10	10
7	14	14	14	14	14	12	12	12	10	10	10	8
8	14	14	14	14	12	12	12	10	10	10	8	8
9	14	14	14	12	12	12	10	10	10	8	8	8
10	14	14	14	12	12	10	10	10	10	8	8	8
12	14	14	12	12	10	10	10	8	8	8	8	6
14	12	12	12	10	10	10	8	8	8	6	6	6
16	12	12	12	10	10	8	8	8	8	6	6	6
18	12	12	10	10	8	8	8	8	8	8	8	5
20	12	12	10	10	8	8	8	6	6	6	5	5
25	10	10	10	8	8	6	6	6	6	5	4	4
30	10	10	8	8	6	6	6	6	5	4	4	3

Wire gauge chart (230V)												
230 Volts	Distance in Feet to Center of Load											
AMPS	20	30	40	50	60	70	80	90	100	120	140	160
5	14	14	14	14	14	14	14	14	14	14	14	14
6	14	14	14	14	14	14	14	14	14	14	14	12
7	14	14	14	14	14	14	14	14	14	14	12	12
8	14	14	14	14	14	14	14	14	14	12	12	12
9	14	14	14	14	14	14	14	14	12	12	12	10
10	14	14	14	14	14	14	14	14	12	12	12	10
12	14	14	14	14	14	14	12	12	12	10	10	10
14	12	12	12	12	12	12	12	12	10	10	10	8
16	12	12	12	12	12	12	12	10	10	10	8	8
18	12	12	12	12	12	12	10	10	10	8	8	8
20	12	12	12	12	10	10	10	10	10	8	8	8
25	10	10	10	10	10	10	10	10	8	8	6	6
30	10	10	10	10	10	10	8	8	8	6	6	6

Electrical Installation & Safety

⚠ DANGER!

- **Never use an extension cord!** An extension cord is any component that adds length to the OEM power cord when connecting the cord to a power source.
- **Never use an adapter plug!** An adapter plug alters the OEM plug configuration when connecting the plug to a power source.
- **Always use the correct outlet.** See "NEMA Plug Configurations" below.

NEMA Plug Configurations

60 HZ USE ONLY!

TRUE uses these types of NEMA plugs shown. If you **DO NOT** have the proper outlet, have a licensed electrician verify and install the correct power source.



International (IEC) Plugs Only

International appliances may be supplied with a power cord that will require installation. Install this cord before connecting the appliance to a power source.

NOTICE > International plug configurations will vary by country and voltage.

Installation

Fully seat the power cord into the appliance receptacle until it locks in position. See fig. 1.

Removal

Depress the red button. See fig. 2.

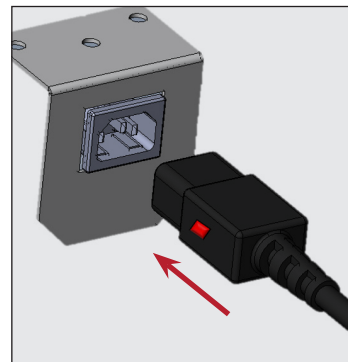


Fig. 1. Fully insert the power cord into the receptacle.



Fig. 2. Push the red button to remove the plug.

About Your Appliance & Installation Requirements (cont.)

Grounded Electrical Duplex with Cover

ⓘ USER ACTION!

Appliances with factory-installed grounded electrical duplex with cover will have two (2) power cords located as described below. **Both power cords require their own separate dedicated circuits.**

- Duplex power cord: Rear or side walls. See fig. 1.
- Appliance power cord: Lower rear. See fig. 2.



Fig. 1. Duplex power cord located on a side wall.

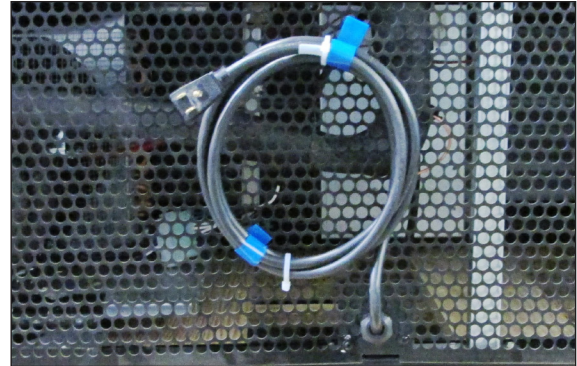


Fig. 2. Appliance power cord located in the lower rear.

Installation & Setup

Installation & Setup

⚠ DANGER!



Risk of electrical shock or burn! Powering off an electronic control or setting temperature controls to the 0 (off) position does not remove power from all components. Unplug the appliance or disconnect power before installation or servicing.

⚠ WARNING!



Only qualified technicians should install and service the appliance. For assistance locating a refrigeration service technician in your area for installation, servicing or maintenance, please visit our Service Company Locator at truemfg.com/support/service-locator/.



The appliance owner is responsible for performing a Personal Protective Equipment (PPE) hazard Assessment and to ensure adequate protection during maintenance and cleaning procedures.

Use appropriate tools, safety equipment, and PPE during installation and servicing.



Sharp edges! Take care when moving, installing, cleaning, servicing and maintaining the appliance to avoid cuts. Be sure to take care when reaching under the appliance or handling metal components. Keep fingers out of pinch point areas, such as the space between appliance doors and surrounding appliances. Take care closing doors with children nearby.



Tip over hazard! Appliance may pose a tipping hazard when uncrating, installing or moving the appliance. Take appropriate safety precautions. Use of tip over restraints may only reduce (not eliminate) the tipping hazard. Never allow children to climb or hang on drawers, doors or shelves.



Crush or cut hazard! Keep clear when uncrating, installing, moving, or servicing the appliance.

Uncrating

ⓘ USER ACTION!



- If the appliance is damaged, note all damage on the delivery receipt, immediately file a claim with the delivery freight carrier, and contact TRUE. **DO NOT install the appliance or put it in service.**
- If the appliance is laid on its back or side, be sure to leave the appliance upright twice the time it has been laying down (up to four (4) hours) before plugging the appliance into a power source. If this time exceeds four (4) hours, let the appliance sit upright for 24 hours before powering the appliance.

ⓘ NOTICE!



Keys for appliances with door locks are located in the warranty packet.

Required Tools

Required tools include (but may not be limited to) the following:

- Adjustable wrench
- Phillips screwdriver

Installation & Setup (cont.)

Procedure

1. Remove the outer packaging (cardboard and bubble wrap or Styrofoam corner and clear plastic). Inspect for concealed damage. See fig. 1. **Again, immediately file a claim with the freight carrier if there is damage.**

NOTICE > **DO NOT** remove the shipping bracket (glass swing doors; fig. 2) or the shipping blocks (glass slide doors; fig. 3) until the appliance is installed in its final location. **Do not discard;** use the bracket/blocks when next moving the appliance.

2. With an adjustable wrench, remove all shipping bolts securing the wood skid to the bottom of the appliance. See fig. 4.

NOTICE > Move the appliance as close as possible to the final location before removing the wooden skid. Some models may require removing the front and/or rear grill/cover to access the shipping bolts.

3. If leveling legs or castors **will not be used**, remove the appliance from the wood skid and set the skid aside.

NOTICE > **DO NOT** lift the appliance by the counter tops, doors, drawers, or grills.

If leveling legs or castors **will be used**, rotate the appliance on the skid (see fig. 5) and see the installation instructions on pg 15.

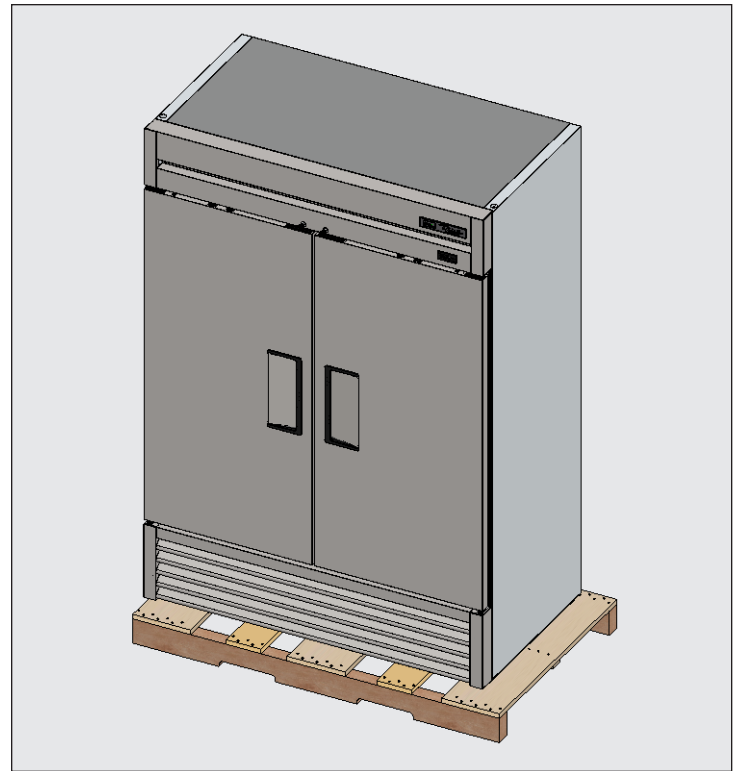


Fig. 1. Remove outer packaging. Inspect for concealed damage.

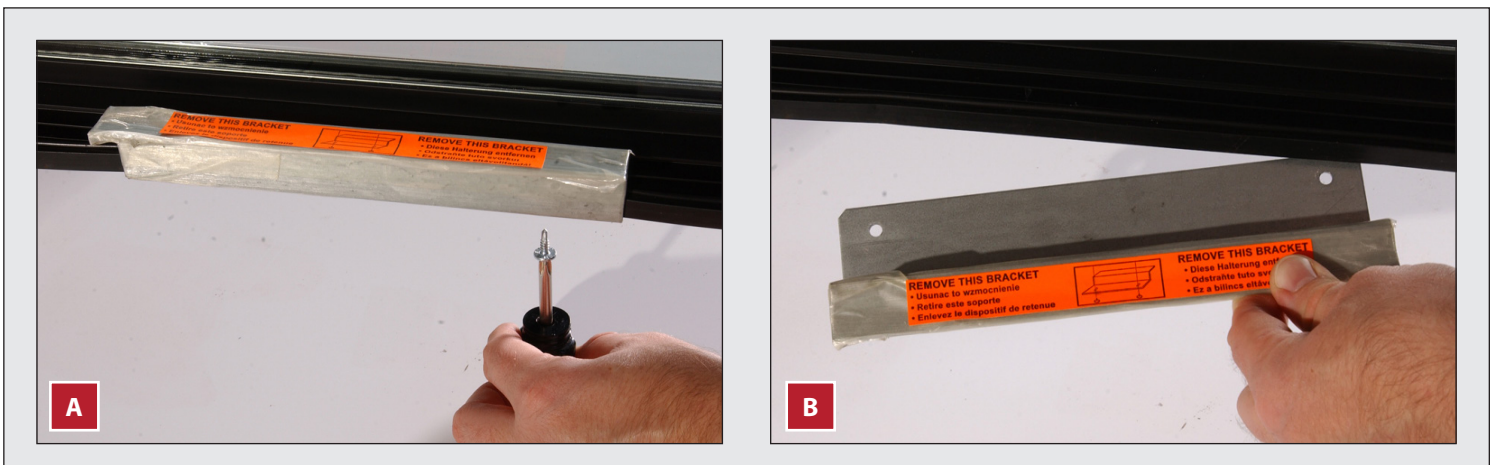


Fig. 2. Remove the glass swing door, if so required. Do not discard.

Installation & Setup (cont.)

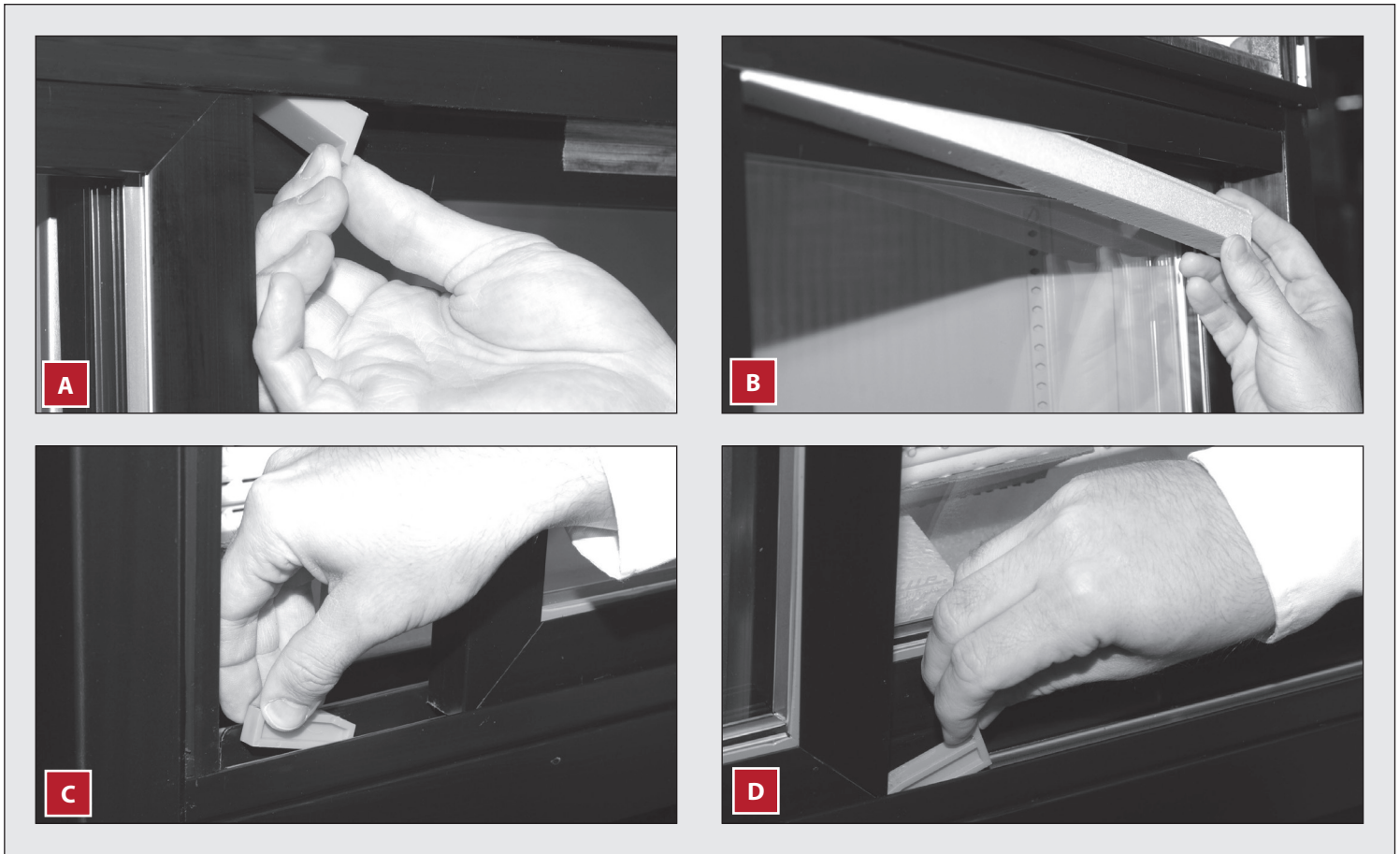


Fig. 3. Glass slide door shipping block locations. Remove if so equipped and do not discard



Fig. 4. Remove the skid bolts.





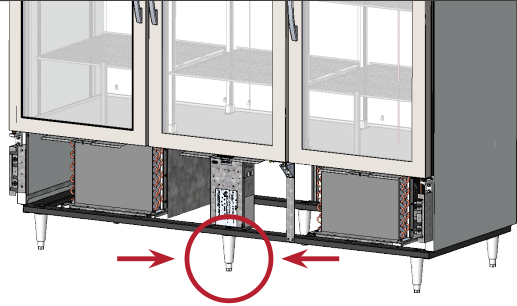

Fig. 5. If leveling legs or castors will be used, rotate the appliance on skid to install.

Installation & Setup (cont.)

Appliance Location

- Ensure that the drain hose or hoses are positioned in the pan.
- Free the plug and cord from inside the lower rear of the appliance (**DO NOT** plug in).
- Place the appliance close enough to the electrical supply so that extension cords are never used.

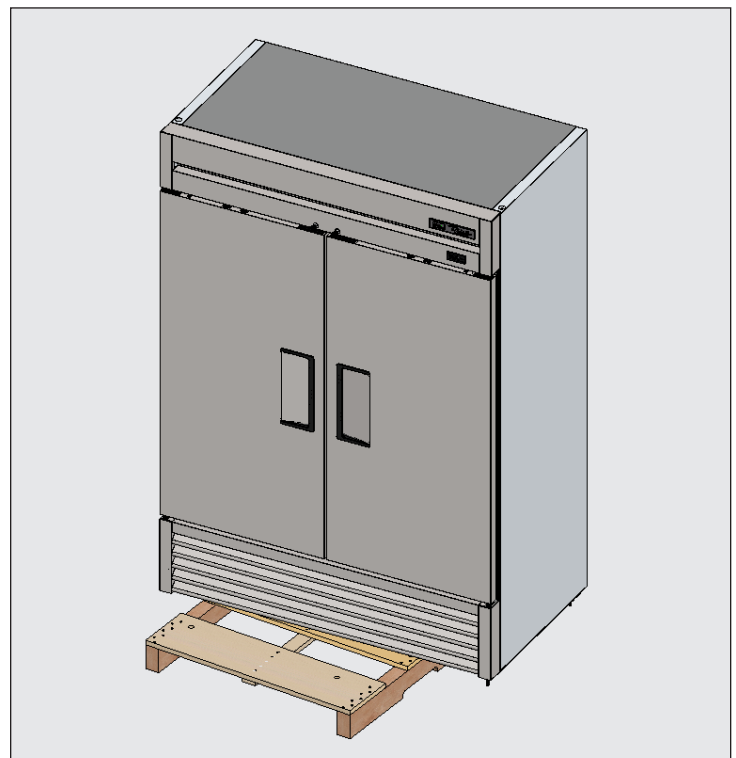Leveling Screw, Leg, or Castor Installation

⚠ WARNING!	
	<p>Tipping hazard! ALWAYS verify center leveling screws fully contact the floor after leveling the appliance.</p> <p>ALWAYS be sure leveling legs or castors (and shims) are snug against the rail assembly or mounting plate.</p> <p>ALWAYS verify center castors or leveling legs fully contact the floor after leveling the appliance. See fig. below.</p>
	
	<p>Sharp edges! Take care when moving, installing, cleaning, servicing and maintaining the appliance to avoid cuts. Be sure to take care when reaching under the appliance or handling metal components.</p>

Leveling screws are provided to assist with leveling the appliance. Adjustable legs provide 6" (152 mm) of clearance under the appliance. Castors provide appliance mobility.

Required Tools

Required tools include (but may not be limited to) the following:

- Adjustable Wrench

Procedure – Leveling Screws

If leveling screws are not factory installed, then, with access to the appliance bottom, install the leveling screws into the threaded holes in the frame rail or appliance bottom. See figs. 1–3.

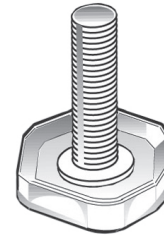


Fig. 1. Leveling screw.

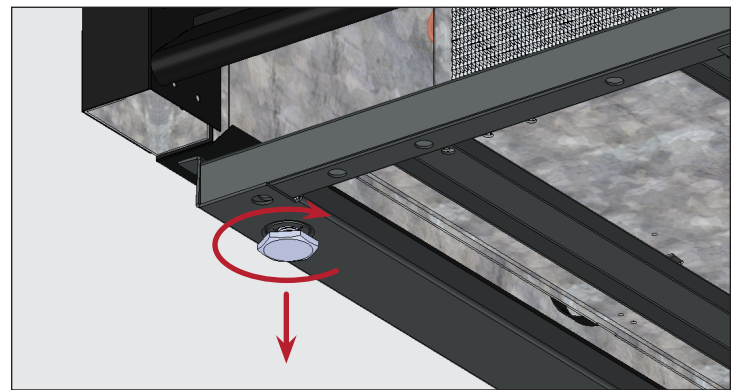


Fig. 2. Turn the leveling screws clockwise to lower the appliance.

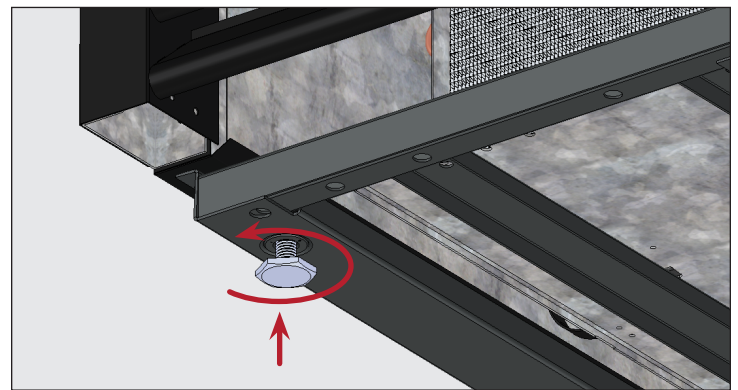


Fig. 3. Turn the leveling screws counterclockwise to raise the appliance.

Installation & Setup (cont.)

Procedure – 6" (152 mm) Leveling Legs

1. Access the appliance bottom.
2. If present, remove the leveling screws.
3. Thread the leveling legs into the frame rail or appliance bottom. See figs. 4 and 5.
4. Verify that the appliance is level. See "Level the Appliance" (pg. 17).
5. If the appliance is not level, gently lift and support the low end of the appliance. Then, with an adjustable wrench, screw the leveling leg bottom stems in or out to level and support the appliance. See fig. 4 and 6.

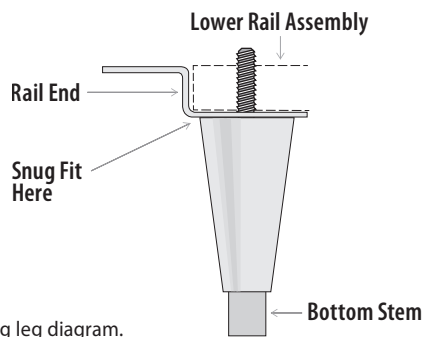


Fig. 4. Leveling leg diagram.

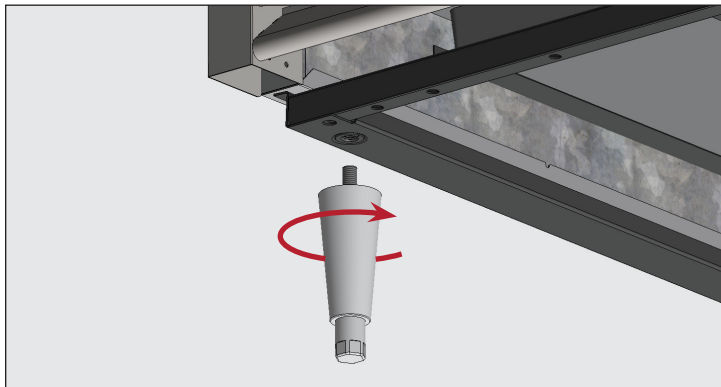


Fig. 5. Screw the leveling legs into the threaded holes.

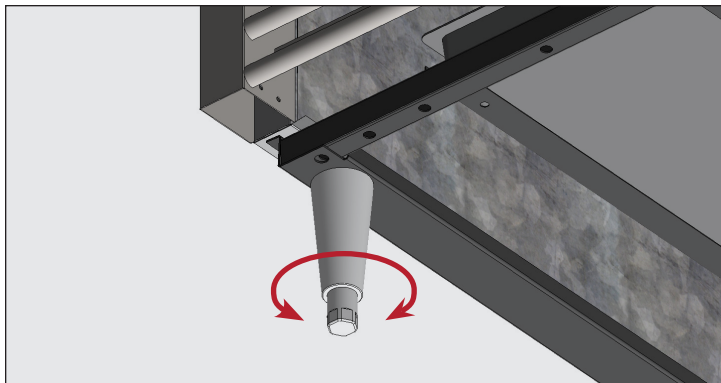


Fig. 6. Turn the bottom stem to level the appliance.

Procedure — Castors

⚠ USER ACTION!



Install castors with brakes in the front.

1. Access the appliance bottom.
2. If present, remove the leveling screws.
3. Thread the castors into the frame rail or appliance bottom. See figs. 7 and 8.
4. Verify the appliance is level. See "Level the Appliance" (pg. 17).
5. On the low end of the appliance, loosen the castor bolt enough to slide provided castor shims between the castor bearing and the bottom rail of the appliance. See fig. 9.
6. Install the desired number of shims. See fig. 9.
 - If more than one shim is used, be sure shim slots are offset. See fig. 10.
 - **DO NOT** use more than four (4) shims on a castor.
 - Be sure each shim touches the castor stem.
7. Tighten and secure the shims and castors.
8. Verify the appliance is level.
9. If the appliance is not level, repeat the steps 3 – 6 until the appliance is level and supported.

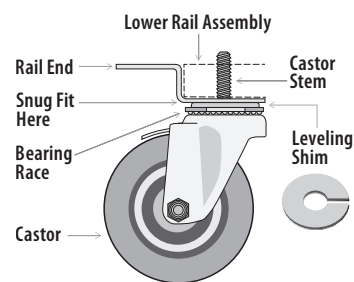


Fig. 7. Castor diagram.

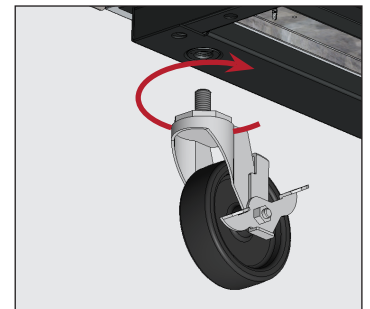


Fig. 8. Screw the castors into the threaded holes.

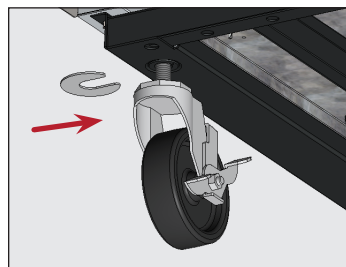


Fig. 9. Insert the castor shim(s), be sure to tighten the castor stem.

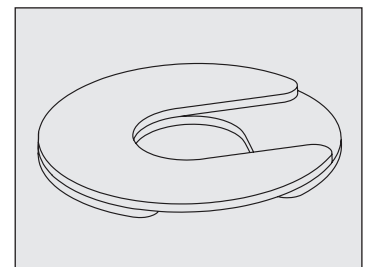



Fig. 10. Install multiple shims with offset slots.

Installation & Setup (cont.)

Level the Appliance

⚠ WARNING!	
	<p>Tipping hazard! ALWAYS verify center leveling screws fully contact the floor after leveling the appliance.</p>

Proper leveling of your TRUE appliance is critical to operating success (for non-mobile appliances). Leveling impacts effective condensate removal and door operation.

See "Leveling Screw, Leg or Castor Installation" (pg. 15) for level adjustment and shimming information.

Procedure

Position the appliance in its final installation location. Then, level the appliance front-to-back and side-to-side. See fig. 1.

1. Position the level on the inside floor of the appliance near the doors (level parallel to appliance front). Level the appliance.
2. Position the level at the inside rear of the appliance (level parallel to appliance back). Level the appliance.
3. Position the level on the left and right inside floor (level parallel to appliance sides). Level the appliance.

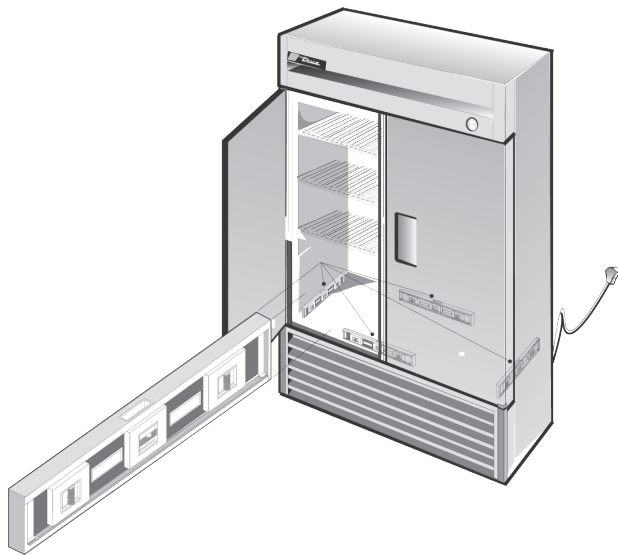



Fig. 1. Measure the level along the perimeter of the interior floor.

Seal the Appliance to the Floor

ⓘ USER ACTION!	
	<p>Asphalt floors are susceptible to chemicals. TRUE recommends installing a layer of tape between the asphalt and the sealant to protect the floor.</p>

The following procedure describes sealing a non-mobile appliance to the floor for NSF standards of proper sanitary operation. Sealing the appliance to the floor prevents liquid spillage on adjacent surfaces of the floor from passing under inaccessible portions of the equipment.

This may not be required for your application.

Procedure

1. Position the appliance in its final installation location. Be sure to leave adequate clearance between the back and sides per "Clearances" (pg. 9) to ensure proper ventilation.
2. Level the appliance front-to-back and side-to-side. See "Level the Appliance" (pg. 17).
3. Draw an outline of the appliance base on the floor.
4. Raise and block the front side of the appliance.
5. Apply a bead of NSF-approved sealant (see list below) to the floor, 1/2" (13 mm) inside the front part of the outline drawn in step 3. The bead of sealant must be heavy enough to seal the entire appliance surface when the appliance is lowered on top of the sealant.
6. Raise and block the rear of the appliance.
7. Apply sealant to the floor on the other three sides, as outlined in step 5.
8. Examine the appliance to ensure that it is sealed to the floor around the entire perimeter.

NSF-Approved Sealants

- 3M #ECU800 Caulk
- 3M #ECU2185 Caulk
- 3M #ECU1055 Bead
- 3M #ECU1202 Bead
- Armstrong Cork – Rubber Caulk
- Products Research Co. #5000 Rubber Caulk
- G.E. Silicone Sealer
- Dow Corning Silicone Sealer

Installation & Setup (cont.)

Shelf Installation

⚠ WARNING!

Overloading, incorrectly installing, or improperly loading shelves can cause shelf failure and negatively affect appliance operation, resulting in appliance damage, product damage, or personal injury.

DO NOT use pliers or any crimping tools when installing shelf clips. Altering shelf clips in any way can lead to shelving instability.



Installation Tips

- Install **all** the shelf clips before installing any shelves.
- Start at the bottom shelf and work your way up.
- Always lay the back of each shelf down on the rear clips before the front.

Procedure

1. Hook the shelf clips into the shelf standards. See fig. 1.
2. Push up on the bottom of the clip. See fig. 2.

NOTICE > You may need to squeeze or twist the bottom of the shelf clip for proper installation. Position all four shelf clips equal in distance from the floor for flat shelves.
3. Ensure the shelf clip is not loose or able to wiggle out of the shelf standard. See figs. 3 and 4.
4. Place the shelves on the shelf clips with the cross support bars facing down.

NOTICE > Be sure all shelf corners are properly seated.

Shelf Adjustment

Shelving is adjustable for customer application. This appliance meets the IEC Shelf Weight Capacity of 47lb/ft² (230kg/m²).

True Track Organizers

True Track Organizers provide a neat and consistent product presentation.



Fig. 1. Installing top tab of shelf clip.



Fig. 2. The bottom tab of the shelf clip will fit tightly.



Fig. 3. You may need to squeeze or twist the bottom of the shelf clip to install.



Fig. 4. Installed shelf clip.

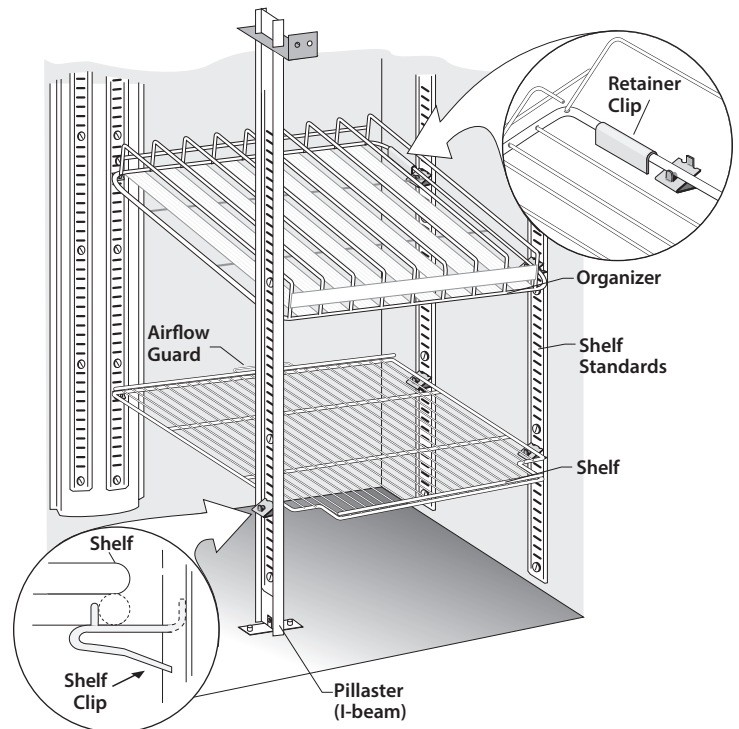


Fig. 5. Close up of a shelf positioned on shelf clips and the retainer clips.

Installation & Setup (cont.)

Chromatography Accessories

Electronic Control (Factory Install Only)

Digital programmable temperature control allows temperature changes in small increments. See fig. 1.

Grounded Electrical Duplex with Cover (Factory Install Only)

- Provides power for stirring devices or other testing equipment inside the appliance.
- Powered with a separate power cord that requires its own dedicated circuit.
- May be mounted on the back, left- or right-inner appliance wall.

Access Port (Factory Install Only)

Access hole in appliance with cover to run lines, cords or move beakers, etc. May be mounted on the back, left- or right-inner appliance wall. See fig. 3.

Chromatography Mast (Factory or Field Install)

Vertical pole inside the appliance to clamp beakers or other equipment to. See fig. 4.



Fig. 1. Electronic temperature control with digital display.



Fig. 2. Grounded electrical duplex with cover.



Fig. 3. Access port.

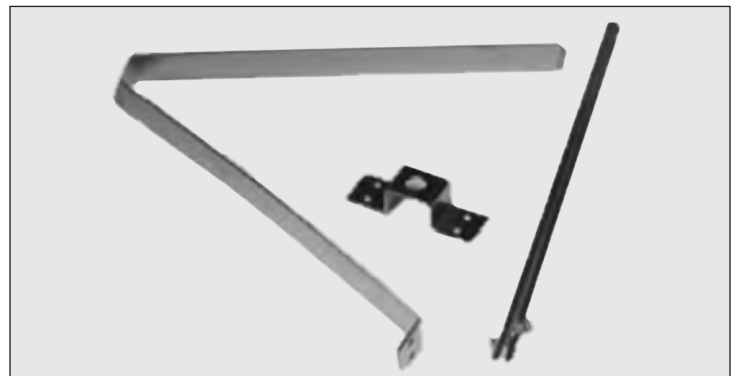








Fig. 4. Chromatography mast components.

Appliance Operation

Appliance Operation

⚠ WARNING – SPOILAGE!	
	<p>CAUTION! The owner is solely responsible for ensuring safe holding temperature levels for perishable items. Failure to do so may result in unsafe product temperatures.</p> <p>Loss or spoilage of products in your appliance is not covered by warranty. In addition to following recommended installation procedures, run the appliance for 24 hours prior to usage to verify operation.</p> <div style="text-align: right;">  </div>



Startup


⚠ DANGER!	
  	<p>Risk of electric shock or fire!</p> <p>Before plugging in the appliance, be sure to inspect the main power cord and plug for damage. Immediately have a qualified technician replace damaged OEM power cords with OEM components.</p>
ⓘ USER ACTION!	
	<p>Before loading product, run your TRUE appliance empty for 24 hours to verify proper operation. Remember, our factory warranty DOES NOT cover product loss!</p>

- The compressor is ready to operate when the appliance is purchased. All you need to do is plug in the appliance.
- Good air flow inside your TRUE appliance is critical. Take care to prevent product from pressing against the sides or back wall and coming within 4" (101.6 mm) of the evaporator housing. Refrigerated air off the evaporator coil must circulate throughout the appliance for even product temperatures.
- Excessive tampering with the control could lead to service difficulties. If replacing the temperature control is ever needed, be sure to order the replacement from your TRUE dealer or recommended service agent.
- All covers and access panels must be in place and properly secured before operating this appliance.

Appliance Operation (cont.)

Temperature Control & Light Switch Location


The light symbol  shows the approximate location of the light switch .

The electronic temperature control can act as a light switch. To control the light, press the up arrow  **M**.



Electronic Temperature Control with Digital Display

On the front top panel.

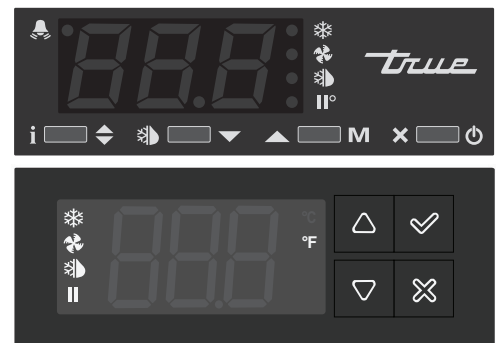

Light Switch on Glass Door Models

Inside the top right ceiling.




Mechanical or Electronic Temperature Control without Digital Display

Inside the top right ceiling or back wall.




Appliance Operation (cont.)

Temperature Control & Light Switch Location (cont.)

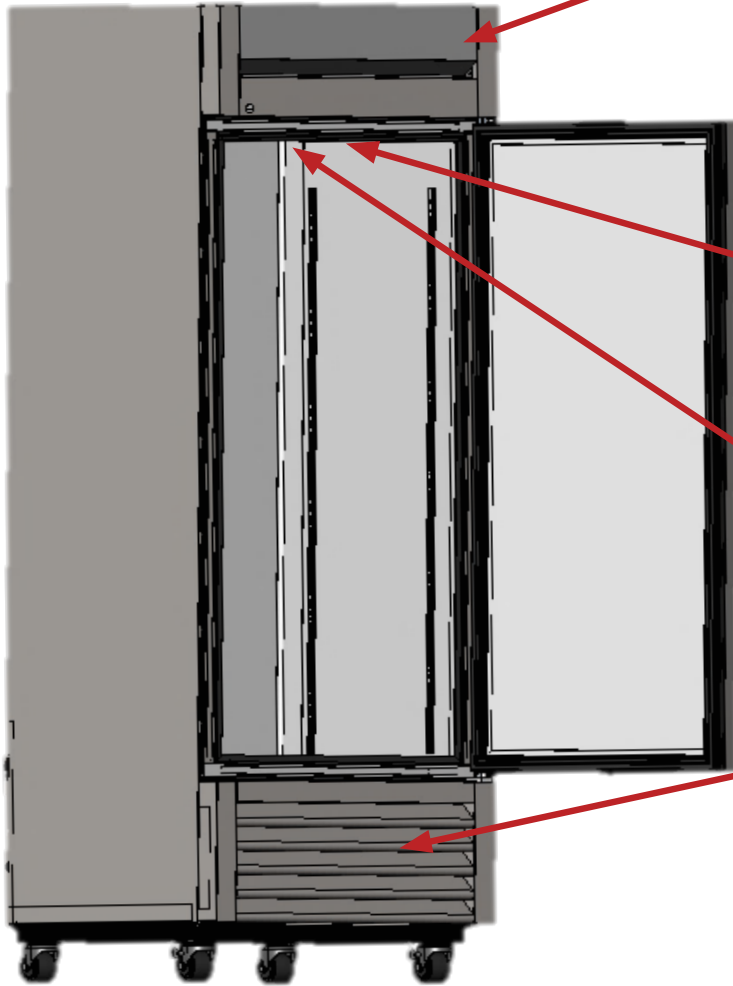


Electronic Temperature Control with Digital Display

On top front of top panel.

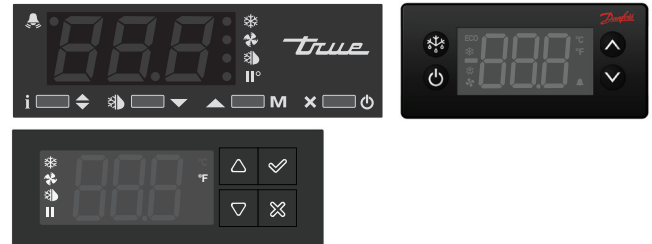


Light Switch on Glass Door Models

Inside the top ceiling.

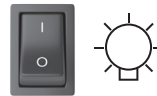

Mechanical Temperature Control

Inside the top ceiling.

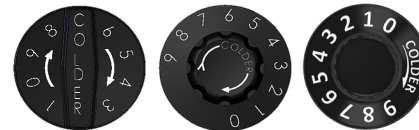

Electronic Temperature Control with Digital Display

Behind the front bottom louvered grill.



Appliance Operation (cont.)

General Sequence of Operation — Refrigerators & Freezers

! USER ACTION!



For more information regarding an appliance's temperature control adjustment or general sequence of operation, please see our Temperature Control Adjustment—Sequence of Operation Manual in our resource library at truemfg.com/support/manuals/#panel4 or scan the QR code.

When the appliance is plugged in...

1. The interior lights illuminate on glass door models (if not, see “Light Switch Location” (pg. 21).
2. The electronic temperature control display illuminates (if installed).

When the appliance is in refrigeration mode

1. There may be a short delay before the compressor and/or evaporator fan(s) start. This delay may be determined by time or by temperature. This delay may also be the result of an initial defrost event that will be a minimum of six (6) min.
2. The temperature control/thermostat may cycle the compressor and evaporator fan(s) on and off together.
3. The temperature control cycles the compressor by sensing either an evaporator coil temperature or air temperature, not a product temperature.

a. Mechanical Temperature Control or Electronic Temperature Control with a Knob

- i. The temperature control cycles the compressor on and off as determined by the cut-in and cut-out temperatures.
 1. #9 is the coldest position
 2. #1 is the warmest position
 3. #0 or Off is the compressor off position

b. Electronic Temperature Control with a Display

- i. The temperature control cycles the compressor on and off as determined by the set point and differential temperatures.
 1. The set point is the adjustable preprogrammed temperature designed to be either the cut-out temperature or it will match the average appliance holding temperature (only when used with a high and low differential).
 2. The differential temperature(s) is the non-adjustable preprogrammed temperature used to determine when the compressor turns on and off (only when the set point is the average holding temperature).
4. An analog thermometer, a digital thermometer, or an electronic control display may reflect the refrigeration cycle swings of up and down temperatures, not a product temperature. **The most accurate method to determine an appliance's operation is to verify the product temperature.**
5. There may be times during refrigeration mode or defrost mode that the condenser fan motor will reverse to blow dirt off the condenser coil.

Appliance Operation (cont.)

General Sequence of Operation — Refrigerators & Freezers (cont.)

When the appliance is in defrost mode...

1. Every appliance requires a defrost event to ensure the evaporator coil remains clear of frost and ice buildup.
2. Defrost is initiated either by the temperature control or a defrost timer.

a. Mechanical Temperature Control

- i. The temperature control cycles the compressor on and off as determined by the cut-in and cut-out temperatures.
 1. During this time, only the evaporator fan runs.
- ii. A freezer with a mechanical temperature control will defrost by time initiation as determined by a defrost timer.
 1. During this time, only the defrost coil heater and drain tube heater are energized.
 2. Defrost terminates when a specific evaporator coil temperature is reached or by a time duration.
- iii. Models with an analog or digital thermometer may show higher-than-normal temperatures during defrost.

b. Electronic Temperature Control

- i. The temperature control is preprogrammed to initiate defrost by a time interval but may also be initiated by temperature demand.
 1. During defrost, a refrigerator turns off the compressor to use the evaporator fans to clear the evaporator coil.
 2. During defrost, a freezer turns off the compressor and evaporator fan to use electric heater to clear the evaporator coil.
- ii. The temperature with a digital display (if installed) shows dEF during defrost.
- iii. Models with an analog or digital thermometer may show higher than normal temperatures during defrost.
- iv. After defrost, there is a display delay until temperature is shown.

NOTICE > The display may have a short delay before showing a temperature after defrost has expired and may show dEF during a refrigeration cycle.

When the appliance sounds an audible and visual alarm...

1. Please reference the appliance's specific temperature control information in Temperature Control Adjustment — Cabinet Sequence of Operation Manual for any alarm codes.

Appliance Operation (cont.)

Temperature Control – **WITHOUT** a Digital Display

See examples below (fig. 1) of temperature control knobs you may find in your appliance.

Turn the temperature control knob clockwise to lower the temperature setting and counterclockwise to raise the temperature setting. One (1) is the warmest setting; Nine (9) is the coldest. See fig. 2.

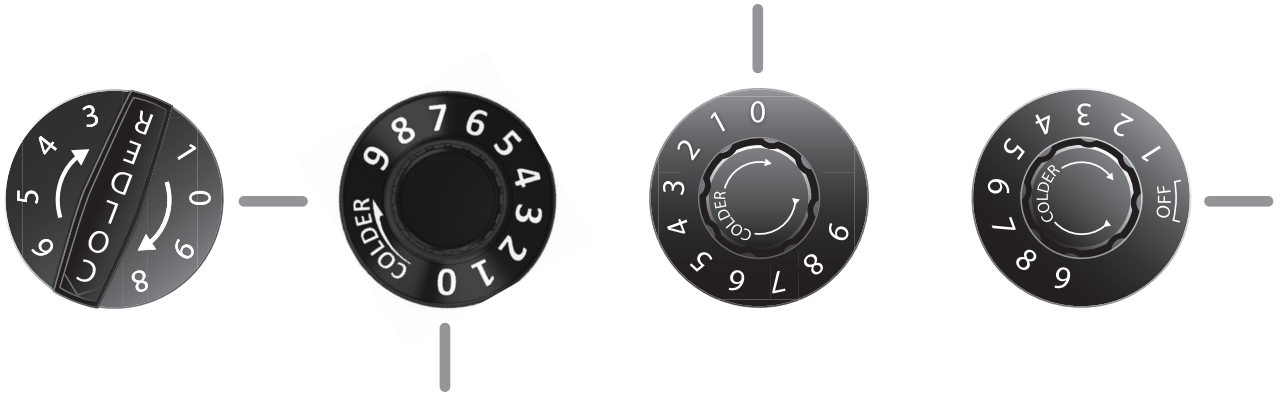


Fig. 1. Temperature control knobs in off position (images not to scale).

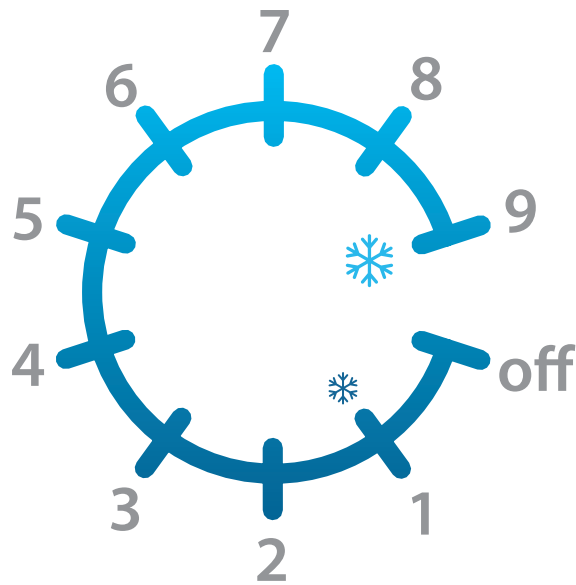


Fig. 2. Temperature control knob settings. 1 is the warmest setting; 9 is the coldest.

Electronic Temperature Control

Electronic Temperature Control

L A E CONTROL LEGEND



- Compressor Running
- Evaporator Fan Running
- Cabinet in Defrost
- Activation of 2nd Parameter Set
- Alarm
- Info/Set Point
- Manual Defrost/Down Arrow
- Manual Activation/Up Arrow
- Stand-By/Cancel

D I S P L A Y C O D E S

Display	Definition	Display	Definition
dEF	Defrost in progress	hi	Cabinet high temperature alarm
oFF	Controller in stand-by	Lo	Cabinet low temperature alarm
do	Door open alarm	E1	Probe t1 failure
t1	Instant probe 1 temperature	E2	Probe t2 failure
t2	Instant probe 2 temperature	E3	Probe t3 failure
t3	Instant probe 3 temperature	thi	Maximum probe 1 temperature recorded
min	Minute of the real time clock	tLo	Minimum probe 1 temperature recorded
hrS	Hours of the real time clock	Loc	Keypad state lock

Electronic Temperature Control (cont.)

LA E CONTROL LEGEND



	Compressor Running		Info/Set Point		Manual Activation/Up Arrow
	Evaporator Fan Running		Manual Defrost/Down Arrow		Stand-By/Cancel
	Appliance in Defrost				
	Activation of 2nd Parameter Set				
	Alarm				

Lock / Unlock the Control

1. Press info to show t1.



2. Press the down arrow until the display shows **Loc**.



3. Press and hold info to show the current lock status.
DO NOT RELEASE THE BUTTON!



YES = Locked



no = Unlocked

4. Press the arrows to change the lock setting.



5. Release all buttons and wait for the control to display the cabinet temperature.

Turn Off / On the Control

⚠ DANGER!

DANGER! Risk of electric shock or burn!
Powering off an electronic does not remove power from all components. Unplug the appliance or disconnect power before installation or servicing.

Turning off the control deactivates all electronic components connected to the control. The lights will remain powered.

Turn Off

Press and hold cancel until the display shows **off**.



Turn On

Press and hold cancel until the display shows **on**.



Electronic Temperature Control (cont.)

LAE CONTROL LEGEND



	Compressor Running		Info/Set Point		Manual Activation/ Up Arrow
	Evaporator Fan Running		Manual Defrost/ Down Arrow		Stand-By/Cancel
	Appliance in Defrost				
	Activation of 2nd Parameter Set				
	Alarm				

Change the Set Point

NOTICE!

The electronic control CANNOT convert a freezer to a refrigerator or vice versa.

Changing the set point adjusts the cabinet temperature to keep optimal product temperature.

1. Press and hold info until the display shows the current set point. **DO NOT RELEASE THE BUTTON!**



2. While holding info, press the up or down arrows to adjust the setting.



3. Release all buttons and wait for the control to display the current cabinet temperature.

Initiate Manual Defrost

A manual defrost clears accumulated frost and ice from the evaporator coil.

The method to initiate a manual defrost is determined by the defrost mode parameter **dFM** setting.

Regular Time Defrost (tiM)

Press and release Manual Defrost until the display shows **dEF**.




Real Time Clock (rtC)

NOTICE > The defrost will only terminate when a specific preset temperature or time duration has been met.

1. Press and hold Manual Defrost until the display shows **dh1**.



2. Press and hold Manual Defrost until the display shows **dEF**.

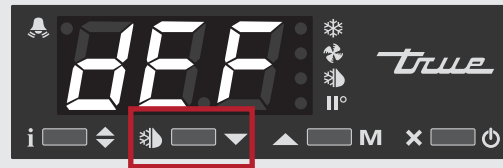

Electronic Temperature Control (cont.)

LAE CONTROL LEGEND



- Compressor Running
- Evaporator Fan Running
- Appliance in Defrost
- Activation of 2nd Parameter Set
- Alarm
- Info/Set Point
- Manual Defrost/Down Arrow
- Manual Activation/Up Arrow
- Stand-By/Cancel

Change System of Measure

The display can show the temperature in either Fahrenheit or Celsius.

1. Press and hold both info and cancel until the display shows **MdL**.



2. Press the down arrow until the display shows **SCL**.



3. Press and hold info until the display shows the current system of measure. **DO NOT RELEASE THE BUTTON!**
 • °F: Fahrenheit • 1°C: Celsius



Fahrenheit



Celsius

4. While holding info, press the up or down arrow to change the system of measure.



5. Release all buttons and wait for the control to display the current cabinet temperature.

Electronic Temperature Control (cont.)

LA E CONTROL LEGEND



- Compressor Running
- Evaporator Fan Running
- Appliance in Defrost
- Activation of 2nd Parameter Set
- Alarm
- Info/Set Point
- Manual Defrost/Down Arrow
- Manual Activation/Up Arrow
- Stand-By/Cancel

Display Probe Temperatures

The display shows the temperature probe readings in different locations of the appliance.

NOTICE > The display may show an average appliance temperature and not a specific probe temperature.

1. Press info to show **t1**.



2. Press and hold info to show the current **t1** probe temperature.



3. Release info to show **t2**.



4. Press and hold info to show the current **t2** probe temperature.



5. Release info to show **t3**.

NOTICE > If t3 probe is not activated, **t3** will not appear.



6. Press and hold info to show the current **t3** probe temperature.



Electronic Temperature Control (cont.)

LA E CONTROL LEGEND



	Compressor Running		Info/Set Point		Manual Activation/Up Arrow
	Evaporator Fan Running		Manual Defrost/Down Arrow		Stand-By/Cancel
	Appliance in Defrost				
	Activation of 2nd Parameter Set				
	Alarm				

Adjust Temperature Alarm Ranges

The temperature alarm ranges are adjustable for different customer applications. Please note the values shown are Fahrenheit.

NOTICE > The control will lock after a short period of inactivity. Unlock the display and press the up arrow to scroll to the required parameter. If the control skips past the necessary parameter, press the down arrow to scroll back.

Step 1: Change Alarm Mode (AtM)

1. Press and hold both info and cancel until the display shows **MdL, ScL, or SPL**.



2. While the display shows **MdL, ScL, or SPL**, press and release info until the display shows **AtM**.



3. While the display shows **AtM**, press and hold info to show the current setting. **DO NOT RELEASE THE BUTTON!**



4. While holding info, press the down arrow until the display shows **AbS** (alarm is an absolute number).



5. Release all buttons to show **ALA**. Proceed to step 2.

Step 2: Change Low Alarm Setting (ALA)

1. While the display shows **ALA**, press and hold info to show the current setting.

DO NOT RELEASE THE BUTTON!



2. While holding info, press the up or down arrows to change the current setting.

NOTICE > Be sure the low alarm setting is **at least 1° colder** than the Cut-Out temperature to prevent a false alarm.



3. Release all buttons to show **AhA**. Proceed to step 3.

Electronic Temperature Control (cont.)

LA E CONTROL LEGEND



	Compressor Running		Info/Set Point		Manual Activation/ Up Arrow
	Evaporator Fan Running		Manual Defrost/ Down Arrow		Stand-By/Cancel
	Appliance in Defrost				
	Activation of 2nd Parameter Set				
	Alarm				

Step 3: Change High Alarm Setting (AhA)

1. While the display shows **AhA**, press and hold info to show the current setting.

DO NOT RELEASE THE BUTTON!



2. While holding info, press the up or down arrows to change the current setting.

NOTICE > Be sure the high alarm setting is at least 1° warmer than the Cut-In temperature to prevent a false alarm.



3. Release all buttons. Proceed to step 4.

Step 4: Change Alarm Delay (Atd)

1. Press info until the display shows **Atd**.



2. While the display shows **Atd**, press and hold info to show the current setting. **DO NOT RELEASE THE BUTTON!**



3. While holding info, press the up or down arrows to change the current setting.


NOTICE >



- Alarm delay is measured in minutes.
- Be sure the time delay is at least 5 min to prevent a false alarm.





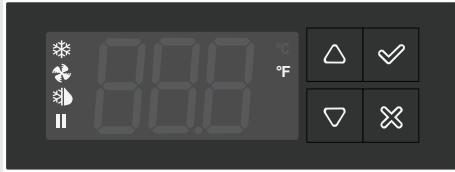



4. Release all buttons and wait for the control to display the current cabinet temperature.

LAE CD25




-  Set point / up arrow
-  Set point / down arrow

-  Confirm / Lights / Info Menu
-  Cancel / Back / Power

-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set

Unlock the Control

1. Press and hold confirm  for 5 seconds. The display will flash **inF** and then show **Loc**.




2. Press confirm to show the current lock status.



YES = Locked



no = Unlocked

3. If **YES**, press the down arrow  to change the setting to **no**.






4. Press confirm.





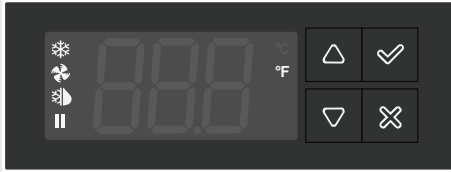



5. The control saves the setting and proceeds to the next item in the info menu. **THE CONTROL WILL LOCK AFTER 60 SEC OF INACTIVITY.**

LAE CD25




-  Set point / up arrow
-  Set point / down arrow

-  Confirm / Lights / Info Menu
-  Cancel / Back / Power

-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set

Turn Off the Control

⚠ DANGER!




Risk of electric shock or burn! Powering off an electronic does not remove power from all components. Unplug the appliance or disconnect power before installation or servicing.

Turning off the control deactivates the refrigeration system and all electronic components connected to the control. The lights will remain powered.

1. Unlock the control. See "Unlock the Control" section. Once unlocked, the display shows **SCL**.

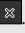

2. Press cancel . The display shows the current appliance temperatures.




3. Press and hold cancel until the display shows **OFF**.



Turn On the Control

1. Unlock the control. See "Unlock the Control" section. Once unlocked, the display shows **SCL**.



2. Press cancel . The display shows **OFF**.







3. Press and hold cancel until the display shows the current appliance temperature.







LAE CD25



-  Set point / up arrow
-  Set point / down arrow

-  Confirm / Lights / Info Menu
-  Cancel / Back / Power


-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set


Change the Set Point



Changing the set point adjusts the appliance operation to keep optimal product temperature.

1. Unlock the control. See "Unlock the Control" section. Once unlocked, the display shows **SCL**.

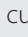

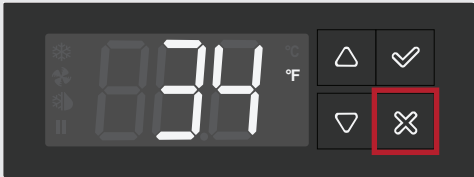2. Press cancel . The display shows the current appliance temperature.



3. Press and hold the up  or down  arrow until the display shows **SP**.



4. With the up or down arrow, adjust the set point to the desired setting.



5. Press confirm , the control saves the setting and the display returns to the current appliance temperature.




Initiate Manual Defrost



A manual defrost clears accumulated frost and ice from the evaporator coil. The defrost will only terminate when a specific preset temperature or duration has been met.

1. Unlock the control. See "Unlock the Control" section. Once unlocked, the display shows **SCL**.

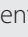

2. Press cancel . The display shows the current appliance temperature.







3. Press and hold both the up  and down  arrow until the display shows **DEF**.







LAE CD25



-  Set point / up arrow
-  Set point / down arrow


-  Confirm / Lights / Info Menu
-  Cancel / Back / Power


-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set

Change System of Measure



1. Unlock the control. See "Unlock the Control". Once unlocked, the display shows **ScL**.

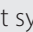

2. Press Confirm  to show the current system of measure.

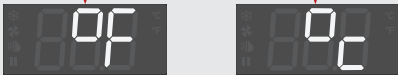

°F = Fahrenheit °C = Celsius

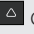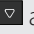3. Press the up  or down  arrow to change the system of measure.




4. With the display showing the desired setting, press confirm.





5. The control saves the setting and proceeds to the next item in the info menu.

Display Probe Temperatures

1. Press and hold confirm  for 5 sec. The display will flash **inF** and then show **Loc**.

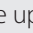

2. Press the up  or down  arrow until the display shows **t1**. This is the thermostat probe.



3. Press confirm. The display shows the current t1 reading.




4. Press confirm. The display shows **t2**. This is the defrost probe.



5. Press confirm. The display shows the current t2 reading.



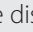



6. Press cancel  twice (x2). The display shows the current appliance temperature.







LAE CD25



-  Set point / up arrow
-  Set point / down arrow


-  Confirm / Lights / Info Menu
-  Cancel / Back / Power

-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set

Temperature Alarms

- The alarm values shown may vary by model. Please see “Adjust Temperature Alarm Ranges” (pg. 38) to adjust the temperature alarms for your application.
- During alarms, the display flashes between the alarm and the current temperature.
- The high temperature alarm has a delay for the unit’s initial startup.
- If an alarm has been silenced but the reason for the alarm persists, the control will sound a 20 sec alert every 60 min until the alarm condition has been resolved as described.

High Temperature Alarm

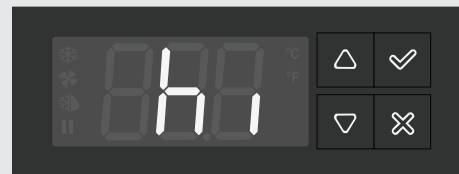

An audio/visual alarm that warns when the internal appliance temperature rises above accepted parameters.


Why it Occurs

If the internal appliance temperature reads $\geq 7^{\circ}\text{C}$ (45°F) the control emits a series of beeps and the display flashes **hi**.

Silencing the Alarm

Press **ANY** button to silence the alarm. The display will continue to flash **hi** until the internal appliance temperature reads $< 7^{\circ}\text{C}$ (45°F).

Low Temperature Alarm

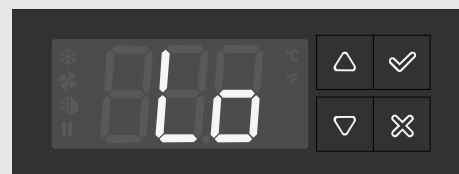

An audio/visual alarm that warns when the internal appliance temperature drops below accepted parameters.


Why it Occurs



If the internal appliance temperature reads $\leq 2^{\circ}\text{C}$ (36°F) the control emits a series of beeps and the display flashes **Lo**.



Silencing the Alarm

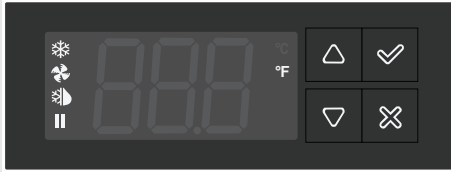



Press **ANY** button to silence the alarm. The display will continue to flash **Lo** until the internal appliance temperature reads $> 2^{\circ}\text{C}$ (36°F).

LAE CD25



-  Set point / up arrow
-  Set point / down arrow

-  Confirm / Lights / Info Menu
-  Cancel / Back / Power

-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set

Door Open Alarm



An audio/visual alarm that warns when the door remains open longer than accepted parameters.

Why it Occurs

If the door remains open for ≥ 60 sec, the control emits a series of beeps and the display flashes **do**.

Silencing the Alarm



Shut the door to stop the alarm. Alternately, press **ANY** button to silence the alarm. The display will continue to flash **do** until the door closes.

Adjust Temperature Alarm Ranges


Step 1: Enter Password

1. Unlock the control. See "Unlock the Control". Once unlocked, the display shows **ScL**.



2. Press the up  or down  arrow until the display shows **PSd**.

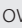


3. Press confirm . The display shows **000**.



4. Press the down arrow until the display shows **989**.




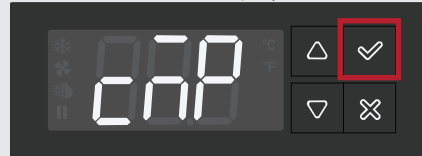

5. Press confirm. The display shows **cMP**.





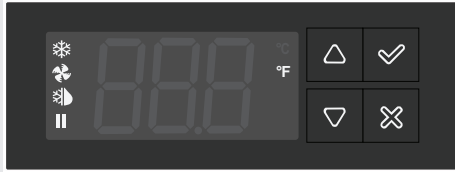



6. Proceed to the next step.

LAE CD25



-  Set point / up arrow
-  Set point / down arrow



-  Confirm / Lights / Info Menu
-  Cancel / Back / Power


-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set


Step 2: Change Alarm Mode (AtM)


1. Once the password is entered, the display shows **cMP**.



2. Press the up  or down  arrows until the display shows parameter group **ALr**.

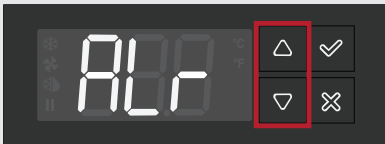

3. Press confirm  to select the parameter group.

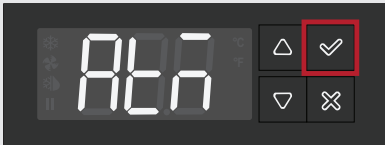


4. Press confirm to select **AtM**. The display shows the current parameter setting.



5. Press the up or down arrow until the display shows **AbS** (alarm is an absolute number).




6. Press confirm to save the setting change. The display shows the next parameter.

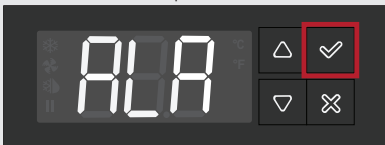


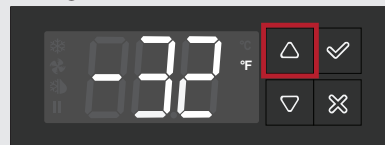
7. Proceed to the next step.

Step 3: Change Low Alarm Setting (ALA)

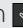
1. Press confirm  to select **ALA**. The display shows the current parameter setting.



2. Press the up  arrow until the display shows the desired setting.



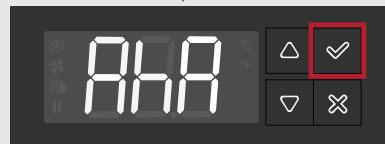

3. Press confirm to enter save the setting change. The display shows the next parameter.









4. Proceed to the next step.

LAE CD25




-  Set point / up arrow
-  Set point / down arrow


-  Confirm / Lights / Info Menu
-  Cancel / Back / Power

-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set

Step 4: Change High Alarm Setting (AhA)


1. Press confirm  to select **AhA**. The display shows the current parameter setting.



2. Press the down arrow  until the display shows the desired setting.




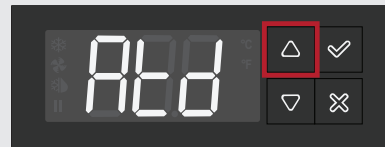
3. Press confirm to enter save the setting change. The display shows the next parameter.

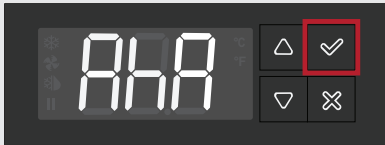


4. Proceed to the next step.


Step 5: Change Temperature Alarm Delay (Atd)


1. Press the up arrow  until the display shows **Atd**.




2. Press confirm  to select **Atd**. The display shows the current parameter settings.




3. Press the down  arrow until the display shows the desired setting.

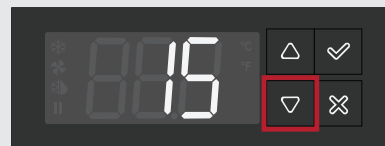


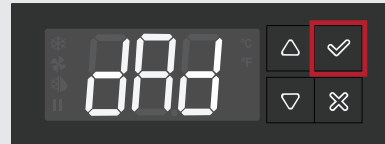

4. Press confirm to enter save the setting change.

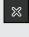



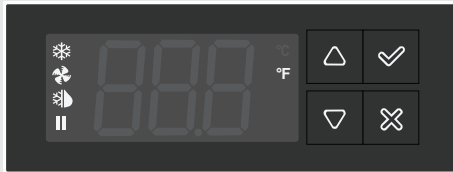



5. Exit the programming.
 - Press cancel  twice (x2)
 - Leave the control inactive for 30 sec

LAE CD25



-  Set point / up arrow
-  Set point / down arrow

-  Confirm / Lights / Info Menu
-  Cancel / Back / Power

-  Compressor Running
-  Evaporator Fan Running
-  Appliance in Defrost
-  Activation of 2nd Parameter Set


High (thi) & Low (tLo) Temperature Log

! USER ACTION!



- The control typically records the **thi** during the initial startup. Reset the **thi** log after 24 hours of operation.
- Reset the log after a high or low temperature alarm.

The high/low temperature log allows users to see the maximum and minimum temperatures record since the log was the last reset.

View the Log

- Press and hold confirm  for 5 sec. The display will flash **inF** and then show **Loc**.



- Press the up  or down  arrow until the display shows **thi** or **tLo**.




- Press confirm. The display shows the logged temperature.





Reset the Log

- Unlock the control. See "Unlock the Control". Once unlocked, the display shows **SCL**.

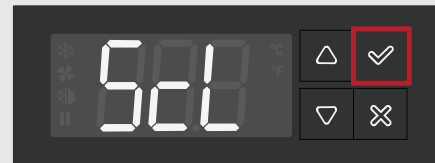


- Press the up  or down  arrow until the display shows **thi** or **tLo**.

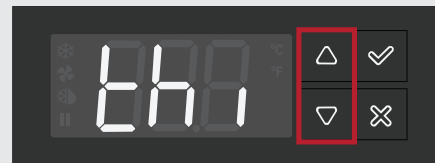

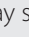- Press confirm . The display shows the logged temperature.



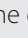





- Press and hold cancel  until the display shows the current internal appliance temperature.






Maintenance & Servicing

Maintenance & Servicing

⚠ DANGER!	
	<p>Risk of electric shock or burn!</p> <ul style="list-style-type: none"> Powering off an electronic control or setting temperature controls to the 0 (off) position DOES NOT remove power from all components. Unplug the appliance or disconnect power before installation or servicing. DO NOT clean appliance with a pressure washer or hose.
	<p>Flammable refrigerant/insulation used! Have a licensed service provider service your appliance to minimize the risk of possible ignition due to incorrect parts or improper service and to ensure the operator's health and safety.</p>

⚠ WARNING!	
	<ul style="list-style-type: none"> Only qualified technicians should install and service the appliance. For assistance locating a refrigeration service technician in your area for installation, servicing or maintenance, please visit our Service Company Locator at truemfg.com/support/service-locator. Turn off and lockout all utilities (gas, electric, water) according to approved practices during maintenance or servicing.
  	<p>The appliance owner is responsible for performing a Personal Protective Equipment (PPE) Hazard Assessment and ensuring adequate protection during maintenance and cleaning procedures.</p> <p>Use appropriate tools, safety equipment, and PPE during installation and servicing.</p>

⚠ WARNING!	
	<p>Sharp Edges!</p> <ul style="list-style-type: none"> Take care when moving, installing, cleaning, servicing, and maintaining the appliance to avoid cuts. Be sure to take care when reaching under the appliance or handling metal components. Stay clear of pinch point areas, such as the space between appliance doors and surrounding cabinetry. Take care closing doors with children nearby.
	<p>Crush or cut hazard! Keep clear when uncrating, installing, moving, or servicing the appliance.</p>
	<p>Slippery Surfaces! Moisture from improper drainage can create slippery surfaces near the appliance. It is your duty to immediately warn your customers of, and dry, the slippery surface. All wet floor areas must be marked with a wet floor sign.</p>

Maintenance & Servicing (cont.)

NOTICE!



- **The appliance owner is responsible** for maintaining the appliance as described in the installation manual. Routine care and maintenance procedures are not covered by TRUE's warranty.
- For additional maintenance instruction, please visit the media center at truemfg.com.
- Any appliance adjustments are to be made **AFTER** the appliance has been verified level and properly supported.

Component Replacement

USER ACTION!



- **Replace components with** original equipment manufacturer (OEM) components such as those shown in fig. 1. OEM parts minimize the risk of possible ignition due to incorrect parts. True is not responsible for defects or damage caused by parts not approved by TRUE. Warranty will be voided for any damage caused by a non-OEM part.
- **Have a licensed service provider** service your appliance to minimize the risk of possible ignition due to incorrect parts or improper service and to ensure the operator's health and safety.



Fig. 1. Example of OEM parts.

Maintenance & Servicing (cont.)


Recommended Maintenance

See recommended maintenance tasks and frequencies below. Some tasks may be required more frequently based on your installation.


Maintenance Tasks	Monthly	Quarterly	Annually
Verify the appliance maintains product temperature.	X	X	X
Inspect power cord for damage; if damaged, replace immediately.	X	X	X
Verify the power cord is fully plugged into the wall receptacle			X
Inspect the overall condition of the appliance and its components (such as castors, doors, and hinges).	X	X	X
Verify operation of all moving parts (such as fan motors, doors and door cords).			X
Check physical condition of all gaskets; verify gaskets seal correctly.		X	X
Inspect any lamps, lamp holder connections, LED modules, and LED module connections.	X	X	X
Check all condenser coils (fronts and backs) for dust and debris; if present, clear the debris.	X	X	X
Check physical condition of all condenser coils and evaporator coils; straighten coil fins as needed.		X	X
Check all evaporator coils for dust and debris; if present, clear the debris.		X	X
Verify the drain line is clear of debris.		X	X


Maintenance & Servicing (cont.)

Condenser Coil Cleaning

⚠ DANGER!	
	<p>Risk of electric shock or burn!</p> <ul style="list-style-type: none"> • Unplug the appliance or disconnect power before installation or servicing. • DO NOT clean appliance with a pressure washer or hose.

⚠ WARNING!	
	<p>The appliance owner is responsible for performing a Personal Protective Equipment (PPE) Hazard Assessment and ensuring adequate protection during maintenance and cleaning procedures.</p>
	<p>Use appropriate tools, safety equipment, and PPE during installation and servicing.</p>
	<p>Sharp edges! Coil fins are sharp and metal components can have sharp edges. Take care when moving, installing, cleaning, servicing, and maintaining the appliance to avoid cuts.</p>
	<p>Risk of eye injury! Airborne dust and debris can cause eye injury. Eye protection recommended.</p>

ⓘ USER ACTION!	
	<p>DO NOT place any filter material in front of the condenser coil.</p>

ⓘ NOTICE!	
	<p>The cleaning of the condenser coil is not covered by warranty!</p>

Required Tools

Required tools include (but may not be limited to) the following:

- Gloves
- Eye Protection
- Phillips Screwdriver or Bit Driver
- 3/8" Driver
- Stiff Bristle Brush
- Tank of Compressed Air
- Vacuum Cleaner
- Drill (optional)

Procedure

1. Unplug the unit or disconnect power
2. Remove the front louver grill. See fig. 1-4.
3. With a stiff bristle brush, carefully clean accumulated dirt from the front condenser coil fins. See fig. 5.
4. With dirt removed from the surface of the coil, use a flashlight to verify you can see clearly through the coil. See fig. 6.

If the view is clear, carefully vacuum any dirt around or behind the condensing unit area. Then reinstall the front louver grill, restore power, and verify operation.

If the view is still blocked with dirt, proceed to the next step.
5. Remove the condenser base bolts. See fig. 7.
6. Remove the condenser coil brackets (if applicable).
7. Carefully slide the condensing unit out (tubing connections are flexible).
8. Gently blow compressed air or CO₂ through the coil until the coil is clean.
9. Carefully vacuum any dirt around and behind the condensing unit area.
10. Reinstall the coil brackets (if applicable). Then, carefully slide the compressor assembly back into position and reinstall.
11. Reinstall the front louver grill, restore power to the unit, and verify operation.

Maintenance & Servicing (cont.)

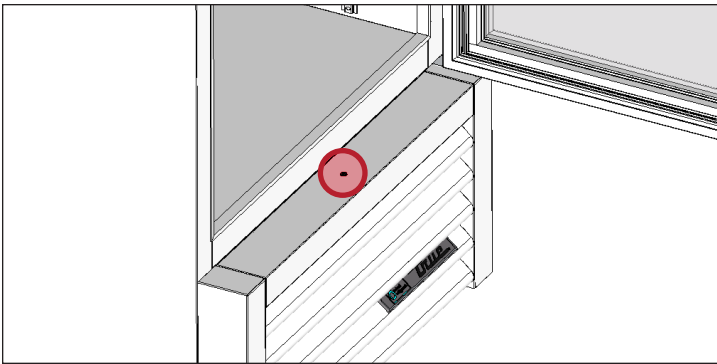


Fig. 1. Swing door: Front louver grill screw location.

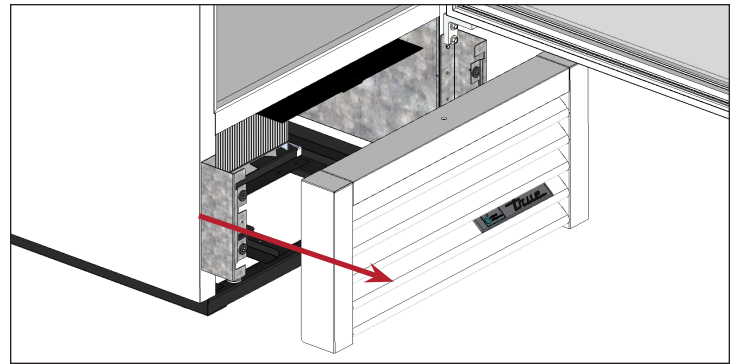


Fig. 2. Swing door: Pull the front louver grill off the appliance.



Fig. 3. Slide Door: Front louver grill screw locations.



Fig. 4. Slide Door: Tilt the top of the grill forward and then lift the grill.

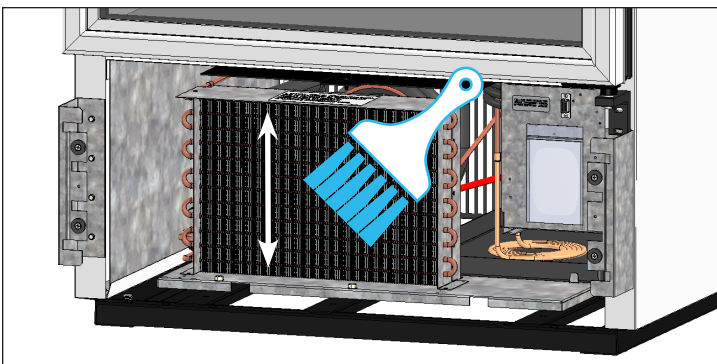


Fig. 5. Never brush across coil fins.

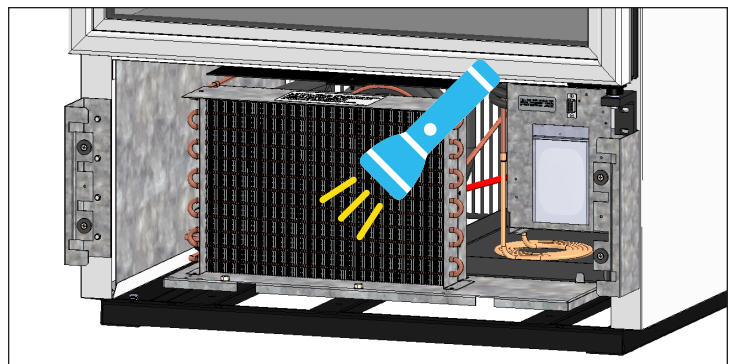


Fig. 6. Verify all blockages have been removed.

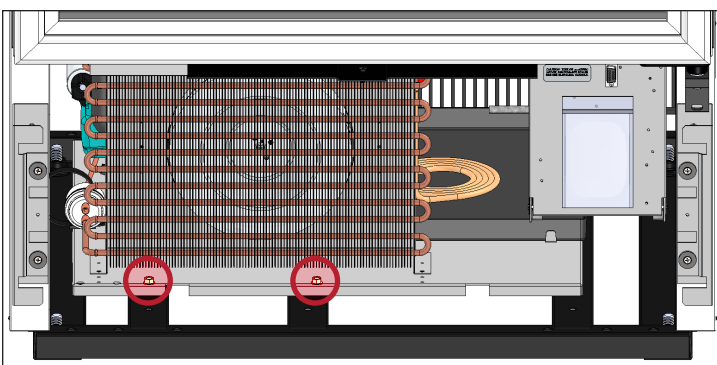





Fig. 7. Condenser base bolt locations.

Maintenance & Servicing (cont.)

General Surface Care & Cleaning

⚠ DANGER!	
	<p>Risk of electric shock or fire!</p> <ul style="list-style-type: none"> • DO NOT clean appliance with a pressure washer or hose.

⚠ WARNING!	
	<p>Slippery Surfaces! Moisture from improper drainage can create slippery surfaces near the appliance. It is your duty to immediately warn your customers of, and dry, the slippery surface. All wet floor areas must be marked with a wet floor sign.</p>
  	<p>The appliance owner is responsible for performing a Personal Protective Equipment (PPE) Hazard Assessment and ensuring adequate protection during maintenance and cleaning procedures.</p> <p>Use appropriate tools, safety equipment, and PPE during installation and servicing.</p>

ⓘ USER ACTION	
	<p>DO NOT USE CITRUS-BASED CLEANERS ON GLASS DOORS.</p>

Gaskets

- Clean gaskets with warm soapy water.
- **DO NOT** use sharp tools or knives to scrape a gasket.
- Avoid full-strength cleaning products.

Glass

- Clean glass with a mild glass cleaner. **DO NOT** use citrus-based cleaners.

Interior

- Clean interior surfaces with a mild solution of baking soda and water to help reduce odor; **DO NOT** use harsh or abrasive cleaners.
- For plastic or powder-coated parts, use warm soapy water to clean **DO NOT** use stainless steel cleaners or similar solvents.

Exterior

- For plastic or powder-coated parts, use warm soapy water to clean **DO NOT** use stainless steel cleaners or similar solvents.
- **DO NOT** clean stainless steel with steel wool or abrasive products. **DO NOT** use detergents or degreasers with chlorides or phosphates. See model-specific manual for more information.

Maintenance & Servicing (cont.)

Stainless Steel Care & Cleaning

! USER ACTION



DO NOT USE ANY STEEL WOOL, ABRASIVE, OR CHLORINE-BASED PRODUCTS TO CLEAN STAINLESS STEEL SURFACES.

Stainless Steel Opponents

There are three basic things which can break down your stainless steel's passivity layer and allow corrosion to form.

- Scratches from wire brushes, scrapers, steel pads, and other items that can be abrasive to stainless steel's surface.
- Deposits left on your stainless steel can leave spots. You may have hard or soft water depending on what part of the country you live in. Hard water can leave spots. Hard water that is heated can leave deposits if left to sit too long. These deposits can cause the passive layer to break down and rust your stainless steel. All deposits left from food prep or service should be removed as soon as possible.
- Chlorides which are present in table salt, food and water, as well as in household and industrial cleaners. These are the worst type of chlorides to use on stainless steel.

Stainless Steel Cleaning and Restoration

Stainless steel cleaners must be free of phosphates, chlorine, chloride, and ammonia.

True offers environmentally-friendly cleaner and polish through our True Store at store.trueresidential.com/products/stainless-steel-clean-polish-kit.

Custom Painted Appliance and Hardware

For painted doors and other surfaces, use a mild solution of soap and water with a soft microfiber cloth.

Maintenance & Servicing (cont.)

8 Tips to Help Prevent Rust on Stainless Steel

Maintain the Cleanliness of Your Equipment

Avoid build-up of hard stains by cleaning frequently. Use cleaners at the recommended strength (alkaline chlorinated or non-chloride).

Use the Correct Cleaning Tools

Use non-abrasive tools when cleaning your stainless steel products. The stainless steel's passive layer will not be harmed by soft cloths and plastic scouring pads.

Clean Along Polishing Lines

Polishing lines ("grain") are visible on some stainless steels. Always scrub parallel to polishing lines when visible. Use a plastic scouring pad or soft cloth when you cannot see the grain.

Use Alkaline, Alkaline-Chlorinated or Non-Chloride Cleaners

While many traditional cleaners are loaded with chlorides, the industry is providing an ever increasing choice of non-chloride cleaners. If you are not sure of your cleaner's chloride content, contact your cleaner supplier. If they tell you that your present cleaner contains chlorides, ask if they have an alternative. Avoid cleaners containing quaternary salts, as they can attack stainless steel, causing pitting and rusting.

Rinse

When using chlorinated cleaners, you must rinse and wipe dry immediately. It is better to wipe standing cleaning agents and water as soon as possible. Allow the stainless steel equipment to air dry. Oxygen helps maintain the passivity film on stainless steel.

Never Use Hydrochloric Acid (Muriatic Acid) on Stainless Steel

Even diluted, hydrochloric acid can cause corrosion, pitting and stress corrosion cracking of stainless steel.

Water Treatment

To reduce deposits, soften hard water when possible. Installation of certain filters can remove corrosive and distasteful elements. Salts in a properly maintained water softener can also be to your advantage. Contact a treatment specialist if you are not sure of the proper water treatment.

Regularly Restore & Passivate Stainless Steel

Stainless steel gets its stainless properties from the protective chromium oxides on its surface. If these oxides are removed by scouring, or by reaction with harmful chemicals, then the iron in the steel is exposed and can begin to oxidize, or rust. Passivation is a chemical process that removes free iron and other contaminants from the surface of stainless steel, allowing the protective chromium oxides to re-form.

Maintenance & Servicing (cont.)

Slide Door Removal

⚠ WARNING!



Crush or cut hazard! Doors are heavy. Be prepared to support the door's weight. Do not lift or support the door by the handle.

🛠 USER ACTION



- **DO NOT** use the side latch before removing the slide door. The tension on the door cord is needed to execute these operation instructions.
- Doors **CANNOT** be removed unless they are positioned as stated in these instructions.

1. Position the doors as described below:
 - **Two-Door Appliances** Slide the front door so it is centered on the appliance. See figs. 1 and 2.
 - **Three-Door Appliances** Center the middle door on the left edge of the right door. See fig. 3.
 2. After centering, lift the door and tilt the top of the door back until the rollers are out of the top channel. Then, swing the bottom of the door out of the bottom channel. See fig. 4.
- Two-Door appliances skip to step 5**
3. Slide the right door to the left and align the left edge with the beginning of the TRUE Logo at the top of the door frame (see fig. 5). Then lift the door out of the track as described in step 2 (see fig. 4).

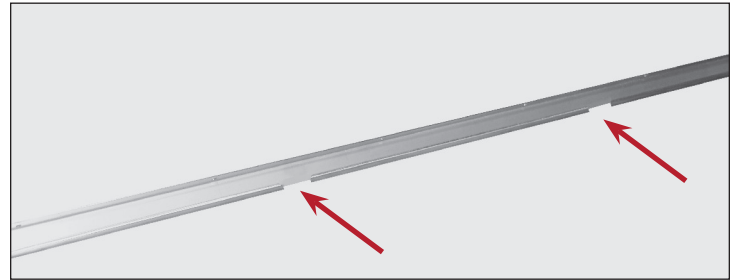


Fig. 1. Door channel openings (two-door appliances ONLY).

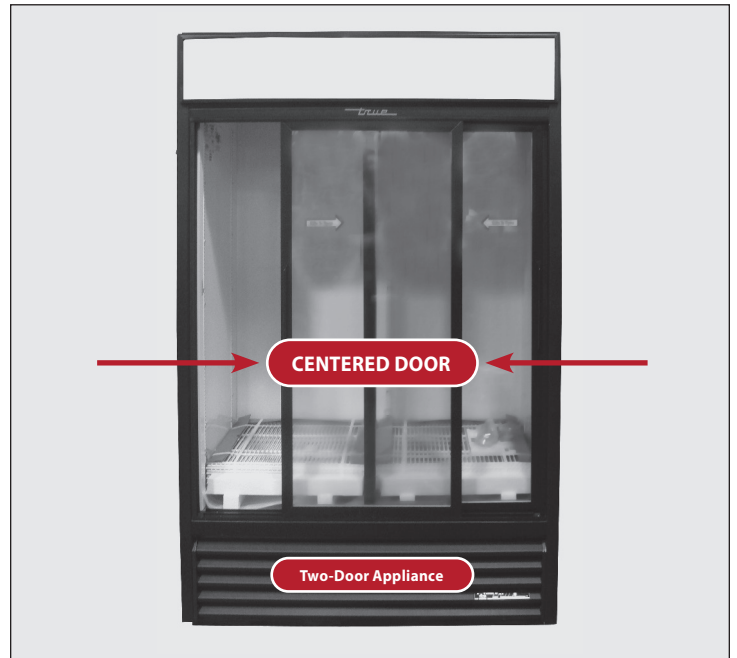


Fig. 2. Position the front door at the center of the appliance.



Fig. 3. Center the middle door on the right door's left edge.



Fig. 4. Carefully remove the door.

Maintenance & Servicing (cont.)

Slide Door Removal (cont.)

1. Slide the left door to the right and align the right edge with the end of the TRUE Logo at the top of the door frame (see fig. 6). Then lift door out of track same way as fig. 4.
2. Remove the door cord from the roller bracket. See figs. 7 and 8.
 - NOTICE >** The door cord will either be nylon cord or metal cable. The black plastic tab holding the door cord slides out the back.
3. Let the door cord slowly retract back into the door side channel.
4. When reinstalling the door, be sure the door cord grommet attaches to roller slot closest to the pulley. See fig. 9.



Fig. 5. Align the right door's left edge with the TRUE logo's left edge.



Fig. 6. Align the left door's right edge with the TRUE logo's right edge.

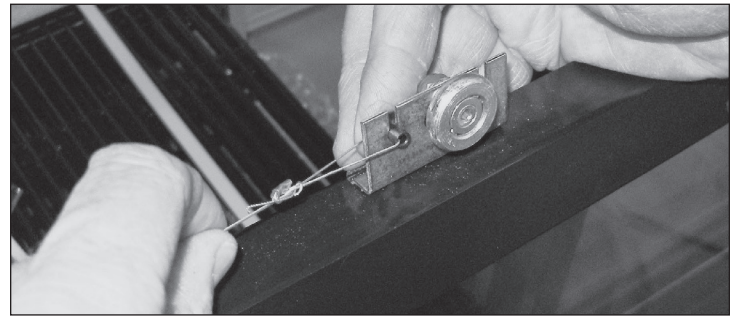


Fig. 7. Remove the door cord from the roller bracket.



Fig. 8. Slide the black tab holding the door cord out the roller bracket's back.

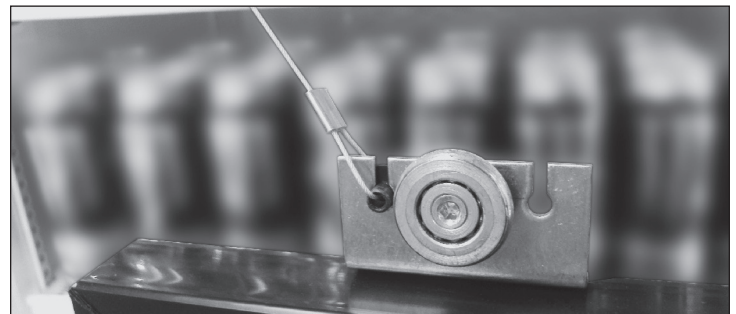


Fig. 9. Be sure the door cord is in the roller slot closest to the pulley.

Maintenance & Servicing (cont.)

Slide Door Adjustment

Only adjust the appliances after it has been installed in its final location and correctly leveled.

Procedure

1. Completely close the slide doors and check for openings. If there are any gaps/openings between the closed doors and appliance, the doors will need to be adjusted.
2. With a 7/16" wrench or adjustable wrench and 1/8" Allen wrench, loosen the roller and move it along the slotted hole. After the adjustment has been made tighten the roller into place. See fig. 1.

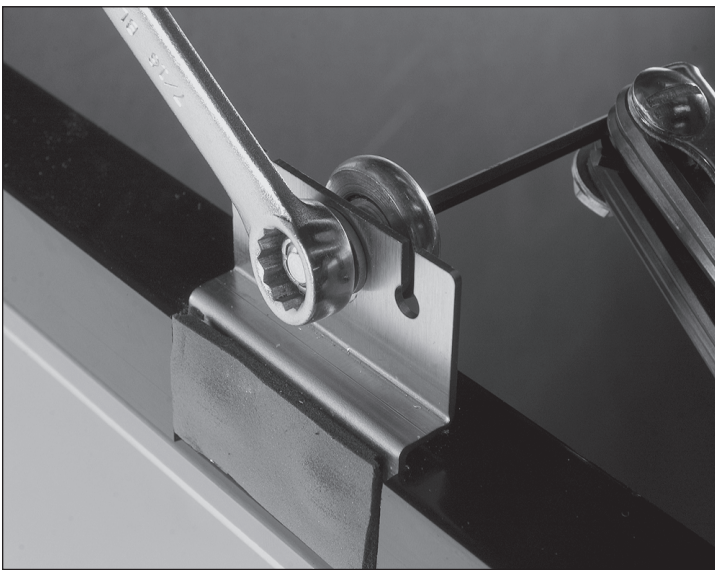


Fig. 1. Adjust the roller.

Slide Door Hold-Open Feature

The hold-open feature keeps the door in the open position.

Procedure

1. Slide the door open.
2. Latch the door in the open position from the back side of door (notch in track). See figs. 1 and 2.

NOTICE > Figs. 1 and 2 show the rear view of the door and track.



Fig. 1. Door latch is in the open position (engaged with notch).



Fig. 2. Door latch is in the closed position (disengaged from notch).

Warranty Information (USA & Canada Only)

Warranty Information

To view and download the
Warranty Information for USA & Canada,
please scan the QR code below.





truemfg.com